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Directions for use | Instrucciones para el uso | Instruction d'emploi | Инструкция по пользованию

Art.-Nr. 0435-9104

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## General safety requirements

- ▶ Read these directions for use carefully before using.
- ▶ Keep these directions for use where they will be accessible for all users.
- ▶ Failure to comply with the directions for use shall render all warranty conditions null and void and may lead to personal injury and material damage.
- ▶ The instrument must only be used by qualified skilled employees.

## Safety instructions

- ▶ Follow the manufacturer's instructions and safety instructions of the medium you are filling.
- ▶ Damage to the respiratory tract and other organs due to off-gassing media. Fill only in well ventilated rooms or under a suitable exhaust hood.
- ▶ The instrument must only be used within the confines of applicable material resistance and for its designated purpose. Any other type of use will affect the warranty/guarantee.
- ▶ The instrument is designed for manual operation.
- ▶ The instrument is not suitable for cleaning in a dishwasher, don't use any sharp cleaners.

## Layout and function

- ▶ The instrument is made of stainless steel V2A (1.4301).
- ▶ The instrument is suitable for safe, fatigue-free filling of liquids. Not suitable for continuous operation.

### NOTE!

The instrument is only suitable for use with Bürkle compact jerrycans 10 litres (see accessories).

- ▶ The device can be used in the temperature range of 0 - 40 °C (briefly up to 60 °C). Do not expose the compact jerrycan to any direct sunlight.
- ▶ The instrument is suitable to fill off fluid and viscosity liquids up to 500 mPas, pay attention to chemical resistance.
- ▶ Constructional changes of the instrument or the accessories aren't allowed.

### NOTE!

- ▶ At maximum filling, take precautions against the liquid running down the jerrycan.

### NOTE!

Abrasion of the Filling support or the compact jerrycan.

- ▶ Before each use do visual control.


## Scope of delivery

1 x	Filling support for compact jerrycan 10 l	
4 x	Plastic feet	⑥

Point out any damage in the packaging to the freight forwarder/supplier immediately.

## Assembly and safe usage

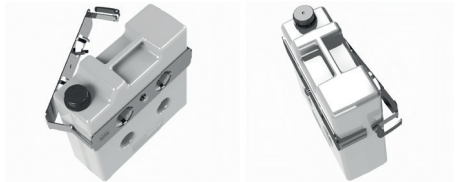1. Attach the holder ① to the closed compact jerrycan 10 l ② and fix it hand-tight with the tightener ③ (see image below).

### CAUTION!

Bodily injury due to squeezing or cutting the fingers.

- ▶ Attach the holder ① carefully.

2. Insert the holder ① with compact jerrycan 10 l in the dispensing stand ④.

3. Tilt opened compact jerrycan 10 l ② at handle ⑤ and fill liquid.

**CAUTION!**

Squeezing and cutting the fingers by uncontrolled swinging of the compact jerrycan ②.

- ▶ Always guide the compact jerrycan ② by the handle ⑤.
- ▶ Keep fingers away from the holder ① and dispensing stand ④ (see pictogram).

**CAUTION!**

Bodily injury due to uncontrolled escaping liquid.

- ▶ Always guide the compact jerrycan ② by the handle ⑤.

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