**Bürkle GmbH**

Rheinauen 5 | D - 79415 Bad Bellingen  
Telefon +49 (0) 76 35 8 27 95-0  
Telefax +49 (0) 76 35 8 27 95-31  
info@buerkle.de | www.buerkle.de

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## General safety requirements

- ▶ Read these directions for use carefully before using.
- ▶ Keep these directions for use where they will be accessible for all users.
- ▶ Failure to comply with the directions for use shall render all warranty conditions null and void and may lead to personal injury and material damage.
- ▶ The instrument must only be used by qualified skilled employees.

## Safety instructions

- ▶ Follow the manufacturer's instructions and safety instructions of the medium you are filling.
- ▶ Damage to the respiratory tract and other organs due to escaping or off-gassing media. Fill only in ventilated rooms or under a suitable exhaust hood. Wear appropriate protective clothing.
- ▶ When working with explosive substances, follow the technical rules for combustible liquids (in accordance with TRBF – (German) Technical Rules on Combustible Liquids) and wear appropriate protective equipment (zone 0 and 1, see annex).

### **WARNING!**

Possible sparking due to electrostatic discharge. Danger of explosion if combustible liquids are used.

- ▶ Thread sealing tape influences the electrical conductivity. The spigot is equipped with a sealing ring. No sealing tape necessary.
  - ▶ Connect the earthing terminal before use (see accessories).
  - ▶ Wear the necessary protective clothing.
- ▶ The instrument must only be used within the confines of applicable material resistance and for its designated purpose. Any other type of use will affect the warranty/guarantee.
- ▶ The instrument is suitable for indoor use and manual operation. Do not expose the device to any direct sunlight.
  - ▶ The instrument is not suitable for cleaning in a dishwasher.

### **CAUTION!**

Escaping liquids due to falling over or tipping of the instrument. Bodily injury due to escaping liquid and damage to the instrument.

- ▶ Make certain the container is positioned with adequate stability.
- ▶ Do not transport with the stopcock screwed on.

## Layout and function

- ▶ The device is made of PA, PTFE and ePTFE.
- ▶ The device is suitable for dispensing non-crystallizing liquids (without solids) from canisters and barrels.
- ▶ The device is suitable for dispensing liquids which require a conductive characteristic, e.g. highly flammable liquids.
- ▶ When filling combustible liquids, always use an antistatic set (see accessories).
- ▶ The device can be used in the temperature range of 0 - 40 °C (briefly up to 60 °C).
- ▶ The device is pressure-resistant up to max. 2 bar at 20 °C.

Point out any damage in the packaging to the freight forwarder/supplier immediately.

### **WARNING!**

Possible electrostatic discharge. Danger of explosion if combustible liquids are used.

- ▶ Connect the earthing terminal before use (see accessories).
- ▶ Wear the necessary protective clothing.

## Assembly and safe usage

### **CAUTION!**

Possible danger of breakage.

- ▶ Apply only normal manual force.
- ▶ Do not lift the container by the stopcock.
- ▶ Do not suspend filling containers from the stopcock.

1. Screw the stop cock with union nut ② onto the container thread (G 3/4" connecting thread), if necessary use thread adapter ③.

### **WARNING!**

Danger of bodily injury due to liquid under high pressure escaping.

- ▶ The pressure in the container must not exceed 2 bar.
- ▶ The discharge reducing nozzle ⑥ is only inserted, when filling with this only pressureless operation is allowed.

### **CAUTION!**

Openings and spaces on the device. Danger of crushing.

- ▶ Be careful for spaces and openings when activating the plug.

2. Open the plug ⑦ (turn anti-clockwise) and dispense liquid.
3. Close the plug ⑦ (turn clockwise).
4. Turn the angle discharge ⑤ up after dispensing to prevent dripping. If necessary, loose the union nut ④, before turning up.

 **WARNING!**

Bodily injury due to escaping liquid

- ▶ Close the stopcock completely after dispensing (set plug to at least „zero“).
- ▶ Keep the stopcock closed when it is not in use.

### Troubleshooting

Problem	To eliminate the error
Thread connection is not sealed	1. Thread is damaged (contact dealer or manufacturer).
Stopcock is dripping	1. Close the plug more tightly. 2. Turn the discharge so it is pointing up.

### Accessories

Type	Specification	Number	Item No.
Antistatic set	to prevent electrostatic charging	1	5602-1000
Compact jerrycan, electroconductive	5 l, with threaded connector 3/4"	1	0435-2055
Compact jerrycan, electroconductive	10 l, with threaded connector 3/4"	1	0435-2105

More accessories can be found at [www.buerkle.de](http://www.buerkle.de)

### Annex

Operational directive in accordance with 1999/92/EC:

Zone 0	a potentially explosive atmosphere is assumed to be continuously present
Zone 1	a potentially explosive atmosphere is assumed to be occasionally present
Zone 2	a potentially explosive atmosphere is not anticipated (and if it does occur, then only rarely and briefly)

Explosion groups:

Group I	for use in mining
Group II	for use in other areas subject to explosion hazard
Group IIA	Maximum permitted gap > 0.9 mm
Group IIB	0.5 mm < maximum permitted gap < 0.9 mm
Group IIC	Maximum permitted gap < 0.5 mm

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