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Directions for use | Instrucciones para el uso | Instruction d'emploi | Инструкция по пользованию

Art.-Nr. 2515-0100

Eyes Fresh Mini-Augendusche

Eyes Fresh mini eyewash station

Mini ducha lavaojos Eyes Fresh

Mini-douche oculaire Eyes Fresh

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General safety requirements

- ▶ Read these directions for use carefully before using.
- ▶ Keep these directions for use where they will be accessible for all users.
- ▶ Failure to comply with the directions for use shall render all warranty conditions null and void.

Safety instructions before using

- ▶ The instrument must only be operated with cold water.
- ▶ Always use and mount the instrument in accordance with its designated use. Any other type of mounting will affect the warranty/guarantee.
- ▶ Use the instrument at room temperature.
- ▶ Do not use the instrument outdoors.

Layout and function

The instrument is freely adjustable. It has dust protection caps ② that drop down automatically during operation. A push button ① on the instrument (arrow) makes it possible to operate the water tap without having to dismount the instrument.

Point out any damage in the packaging to the freight forwarder/supplier immediately.

Assembly

1. If there is a jet controller on the water tap, remove it before assembling the installation.
2. Screw the instrument onto the water tap (standard thread M22x1"). If necessary remove the premounted adapters (IG 3/8"x outer thread 1/2" and outer thread M22xinner thread 1/2").
3. Align the instrument so that the push button is pointing forward.
4. Affix the label with the note on it (included with delivery) next to the wash basin where it is visible.

Operation of the mini eyewash station

CAUTION!

If the stream of water is too strong, it can lead to eye injuries.

- ▶ Adjust the stream of water with the water tap (max. height 30 cm).

1. Open the water tap.
2. Open the eyelids with fingers and flush the eyes with plenty of liquid.
3. If you come in contact with chemicals, consult a physician and follow the directions of the chemical producer.

After the protective dust covers have been moistened, place them back on the nozzles again.

Functional check

1. Open the water tap for 20 s weekly.
2. Close the water tap and place the protective dust covers that had dropped down back on the nozzles.

Maintenance

Check the instrument regularly for impurities (deposits on or in the tube, lime).

1. Unscrew the instrument from the water tap.
2. Remove the protective dust caps.
3. Rinse the instrument with water.
4. If there are heavy impurities, place the instrument in a descaler (follow the directions on the descaler) and then rinse with water.

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