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Directions for use | Instrucciones para el uso | Instruction d'emploi | Инструкция по пользованию

Art.-Nr. 5005-8012

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OTAL® disposable hand pump

Bomba manual desechable OTAL®

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### General safety requirements

- ▶ Read these directions for use carefully before using.
- ▶ Keep these directions for use where they will be accessible for all users.
- ▶ Failure to comply with the directions for use shall render all warranty conditions null and void and may lead to personal injury and material damage.
- ▶ The instrument must only be used by qualified skilled employees.

### Safety instructions before using

- ▶ Follow the manufacturer's instructions and safety instructions of the medium you are filling.
- ▶ Damage to the respiratory tract and other organs due to off-gassing media. Fill only in ventilated rooms or under a suitable exhaust hood.
- ▶ The instrument must only be used within the confines of applicable material resistance and for its designated purpose. Any other type of use will affect the warranty/guarantee.

#### **! WARNING!**

The instrument is not suitable for highly flammable or explosive liquids.

- ▶ The instrument must only be used within the confines of applicable material resistance and for its designated purpose.

#### **NOTE!**

Before using, the instrument has to be cleaned to the extent of your (validated) cleaning process.

- ▶ Measures of manual cleaning have to be carried out in the cleaning/desinfection area and under consideration of occupational health and safety.
- ▶ Pay attention to the material resistance.

### Safety instructions during and after use

- ▶ The instrument generates overpressure. Possible bodily injury due to escaping liquids.
- ▶ The instrument is not suitable for cleaning in a dishwasher.

### Layout and function

- ▶ The disposable riser pipes are made of PE and the universal bungs of PVC.
- ▶ The instrument is suitable for filling thin to low viscosity liquids up to 500 mPas.
- ▶ The instrument is suitable at a temperature range-

from 0 to 40 °C.

#### **! WARNING!**

The system is not tight-closing. Possible gas leakage. Bodily injury due to escaping gases.

- ▶ Always keep valves open when not in use.
- ▶ Ensure adequate ventilation.
- ▶ Wear the necessary protective clothing.

- ▶ The bung fits onto container openings with a diameter of 30 to 70 mm forming an airtight seal.

Ⓐ Ø 30 - 50 mm

Ⓑ Ø 40 - 60 mm

Ⓒ Ø 50 - 70 mm

### Scope of delivery

10 x	Delivery tube, disposable	①
1 x	Universal bungs, three parts	②
1 x	Pump ball	④
1 x	Cutter	③

- ▶ Check delivery for completeness and transport damage.
- ▶ Point out any damage in the packaging to the freight forwarder/supplier immediately.

### Assembly

1. Insert the delivery tube ① into the universal bung ② (opening "Rohr 12 mm") and push it through to the outlet.

#### **NOTE!**

If necessary adjust the universal bung to the opening diameter of the container.

- ▶ When using all three parts (Ⓐ, Ⓑ and Ⓒ) loosen Ⓑ and Ⓒ a bit of part Ⓐ to push the delivery tube through.

2. Determine the conveying tube length on the basis of the container height and shorten the delivery tube ① slightly at an angle to the container bottom using the cutter ③.

**NOTE!**

If necessary, shorten the discharge of the delivery tube ① with the cutter ③ to the desired length.

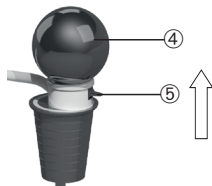
3. Insert the delivery tube ① with the universal bung ② into the container and press it in tightly.
4. Insert the pump ball ④ into the universal bung ② (opening „Pumpe“).

**! WARNING!**

Spraying of liquid under high pressure.

- ▶ Hoses and connections must be faultless.
- ▶ Ensure correct installation.

5. Close the stop-valve ⑤ (push the lever away from the connecting hose).

**Safe filling**

1. Hold the discharge tube above the filling container.
2. Activate the pump ball ④ and generate overpressure. Liquid starts to flow.
3. To complete the filling process open the stop-valve ⑤ (push the lever to the connecting hose) and allow the overpressure to escape.


**Troubleshooting**

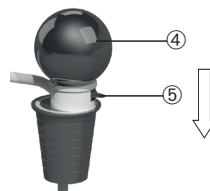Problem	To eliminate the error
The instrument does not generate an overpressure	<ol style="list-style-type: none"> <li>1. Stop-valve is opened.</li> <li>2. The instrument is not connected with the container to form a seal.</li> </ol>
Instrument is not pumping any liquid	<ol style="list-style-type: none"> <li>1. The instrument is not connected with the container to form a seal.</li> <li>2. The riser pipe does not extend into the liquid being filled.</li> <li>3. The delivery tube is not cut at an angle at the height of the tank bottom.</li> </ol>

**Accessories**

Type	Specification	Number	Item No.
Delivery tube PE, disposable	For OTAL® disposable pumps	10	5000-8000

More accessories can be found at [www.buerkle.de](http://www.buerkle.de)

**! WARNING!**

The container is under pressure. Bodily injury due to escaping liquid.

- ▶ Allow excess air to escape completely from the container.
- ▶ After pressure reduction, drain off any remaining liquid of the discharge tube completely.

**NOTE!**

The stop-valve is not suitable for precise dosing.

**Completing filling and dismantling**

1. Pull the pump ball ⑤ off the universal bung ②.
2. Carefully remove the universal bung ② with delivery tube ① from the container.
3. Cut off the delivery tube ① directly below the universal bung ② using the cutter ③.
4. Pull the upper end of the delivery tube ① out of the universal bung ② and dispose it properly with the lower end.

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