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Directions for use | Instrucciones para el uso | Instruction d'emploi | Инструкция по пользованию

Art.-Nr. 5201-1000

Pump-it® Kanisterpumpe

Pump-it® container pump

Bomba para recipiente Pump-it®

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General safety requirements

- ▶ Read these directions for use carefully before using.
- ▶ Keep these directions for use where they will be accessible for all users.
- ▶ Failure to comply with the directions for use shall render all warranty conditions null and void and may lead to personal injury and material damage.
- ▶ The instrument must only be used by qualified skilled employees.

Safety instructions before using

- ▶ Follow the manufacturer's instructions and safety instructions of the medium you are filling.
- ▶ Damage to the respiratory tract and other organs due to off-gassing media. Fill only in ventilated rooms or under a suitable exhaust hood.
- ▶ The instrument must only be used within the confines of applicable material resistance and for its designated purpose. Any other type of use will affect the warranty/guarantee.

WARNING!

The instrument is not suitable for highly flammable or explosive liquids.

- ▶ The instrument must only be used within the confines of applicable material resistance and for its designated purpose.

NOTE!

Before using, the instrument has to be cleaned to the extent of your (validated) cleaning process.

- ▶ Measures of manual cleaning have to be carried out in the cleaning/desinfection area and under consideration of occupational health and safety.
- ▶ Pay attention to the material resistance.

Safety instructions during and after use

- ▶ The instrument generates overpressure. Possible bodily injury due to escaping liquids.
- ▶ The instrument is not suitable for cleaning in a dishwasher.

Layout and function

- ▶ The instrument is made of PP, PE and FKM.
- ▶ The instrument is suitable to fill off highly fluid and low-viscosity liquids.
- ▶ The pump thread is suitable for container threads 64 mm Ø (BSI).
- ▶ The instrument is maintenance-free.

Point out any damage in the packaging to the freight forwarder/supplier immediately.

Assembly

1. Insert the telescopic riser pipe ① into the pump up to the stop.
2. Pull out the telescopic riser pipe completely.
3. Screw the pump onto the container thread. If necessary use a matching thread adapter (see accessories).

Safe filling

1. Hold the discharge tube above the filling container.
2. Activate the manual piston ③ and generate overpressure.
3. To fill off liquid, open the drain valve ② (push the lever downwards).

Completing filling and dismantling

CAUTION!

The container is under pressure. Bodily injury due to escaping liquid.

- ▶ Slowly remove the instrument from the container.

1. Carefully remove the instrument from the container.
2. Rinse the instrument with a pH-neutral fluid.

Troubleshooting

Problem	To eliminate the error
The instrument does not generate an overpressure	1. The instrument is not connected with the container to form a seal.
Instrument is not pumping any liquid	1. The instrument is not connected with the container to form a seal. 2. The riser pipe does not extend into the liquid being filled.

Accessories

Type	Specification	Item No.
Adapterset Pump-it	Consisting of: 2" coarse Mauser, 2" coarse TriSure, 2" BSP, DIN 60, DIN 51	5201-0030

More accessories can be found at www.buerkle.de

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