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Directions for use | Instrucciones para el uso | Instruction d'emploi | Инструкция по пользованию

Art.-Nr. 5305-0110

MiniSampler PTFE

MiniSampler PTFE

MiniSampler PTFE

MiniSampler PTFE

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General safety requirements

- ▶ Read these directions for use carefully before using.
- ▶ Keep these directions for use where they will be accessible for all users.
- ▶ Failure to comply with the directions for use shall render all warranty conditions null and void and may lead to personal injury and material damage.
- ▶ The instrument must only be used by qualified skilled employees.

Safety instructions before using

- ▶ Follow the manufacturer's instructions and safety instructions of the medium you are sampling.
- ▶ When sampling hazardous substances, follow the technical rules on hazardous substances (in accordance with TRGS – (German) Technical Rules on Hazardous Substances) and wear appropriate protective equipment.
- ▶ Damage to the respiratory tract and other organs due to off-gassing media. Use in ventilated rooms or under a suitable exhaust hood.
- ▶ The instrument is not suitable for sampling highly flammable/explosive liquids.
- ▶ The instrument is not suitable for aggressive acids and bases. Vapours could destroy pump.

NOTE!

Before using, the instrument has to be cleaned to the extent of your (validated) cleaning process.

- ▶ Measures of manual cleaning have to be carried out in the cleaning/desinfection area and under consideration of occupational health and safety.
- ▶ Pay attention to the material resistance.

Safety instructions during and after use

CAUTION!

Liquid entering the instrument due to overfilling or tipping. Bodily injury due to escaping liquid and damage to the instrument.

- ▶ Shut down and secure the instrument immediately.
- ▶ Contact the dealer or manufacturer.
- ▶ The instrument must only be used within the confines of applicable material resistance and for its designated purpose. Any other type of use will affect the warranty/guarantee.
- ▶ Do not hold the instrument inclined, lay it down or store it, when the sample box is filled.
- ▶ The instrument is not suitable for cleaning in a dishwasher.

Layout and function

- ▶ The instrument is made of non-conductive material and is suitable for sampling non-combustible liquids from a maximum suction depth of 5 m.
- ▶ The instrument is suitable for sampling liquids from 0 - 60 °C.
- ▶ The instrument is only suitable to take samples and not for continuous operation.
- ▶ Only use the instrument with the recommended sample boxes specified in the accessories.
- ▶ Protect the instrument against effects of weather exposure during storage.
- ▶ The instrument is maintenance-free.

Parts in contact with medium

- ▶ PTFE (suction hose)
- ▶ PFA (suction box)
- ▶ Stainless steel, brass nickel-plated (hose weight)

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Scope of delivery

1 x	MiniSampler Vacuum pump	①
1 x	Sample box, 180 ml (PFA)	②
1 x	Suction hose (PTFE, 5 m)	③
1 x	Hose weight (stainless steel, brass nickel-plated))	④
1 x	Hose cutter	
1 x	Directions for use	
20 x	Seal-it lead bottle seals	
1 x	Transport case	

- ▶ Check delivery for completeness and transport damage.
- ▶ Point out any damage in the packaging to the freight forwarder/supplier immediately.

Assembly

WARNING!

Danger of cutting when using the hose cutter.

- ▶ Keep hands away from the cutting area.

1. Screw the sample box ② onto the instrument finger-tight.
2. Guide the suction hose ③ through the open clamping screw ⑤ about 1 cm into the sample box.
3. Screw on the clamping screw finger-tight and fix the suction hose in place.
4. Place the free end of the suction hose in the liquid you are sampling. Ensure that the suction hose is not kinked.
5. Place the hose weight ④ on the suction hose depending on the location where it will be used and tighten the union nut. The hose weight improves the stability of the suction hose.

NOTE!

Possible loss of the hose weight during use.

- ▶ Screw on the hose weight tightly with the cap nut.
- ▶ Verify that the hose weight is attached correctly before use.

Safe sampling

WARNING!

Bodily injury due to escaping media when overfilling the sample box. Liquids can be aspirated into the piston and can escape at the back of the instrument.

- ▶ Do not overfill the sample box.
- ▶ Wear the appropriate protective clothing.

1. Activate the handle ⑥ and pump liquid into the sample box.
2. Screw off the sample box and seal it.

WARNING!

Bodily injury due to escaping media when screwing off the sample box and pulling the suction hose out of the instrument.

- ▶ Dismount the instrument carefully.
- ▶ Wear the appropriate protective clothing.

After taking the sample

1. Open the clamping screw and remove the suction hose.
2. Rinse the suction hose (together with the hose weight), sample box and instrument with a pH-neutral liquid.

Troubleshooting

Problem	To eliminate the error
Pump is not generating a negative pressure	Repeat assembly steps 1 to 4. Seal or valve faulty (contact dealer or manufacturer).

Accessories

Type	Specification	Number	Item No.
Sample box (PFA)	90 ml	1	6205-0090
Sample box (PFA)	180 ml	1	6205-0180
Suction hose (PTFE, on roll)	Ø 6x8 mm, length 5 m	1	8804-0608
PTFE MiniSampler adapter	for glass thread GL45	1	5305-0209

More accessories can be found at www.buerkle.de

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