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Directions for use | Instrucciones para el uso | Instruction d'emploi | Инструкция по пользованию

Art.-Nr. 5307-1055 / 5307-1085 / 5307-1120 / 5307-2055 / 5307-2085 / 5307-2120

MicroSampler

MicroSampler

MicroSampler

MicroSampler

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General safety requirements

- ▶ Read these directions for use carefully before using.
- ▶ Keep these directions for use where they will be accessible for all users.
- ▶ Failure to comply with the directions for use shall render all warranty conditions null and void and may lead to personal injury and material damage.
- ▶ The instrument must only be used by qualified skilled employees.

Safety instructions before using

- ▶ Follow the manufacturer's instructions and safety instructions of the medium you are sampling.
- ▶ Do not fill the instrument with potentially explosive substances.
- ▶ The instrument must only be used within the confines of applicable material resistance and for its designated purpose. Any other type of use will affect the warranty/guarantee.
- ▶ When sampling hazardous substances, follow the technical rules on hazardous substances (in accordance with TRGS – (German) Technical Rules on Hazardous Substances) and wear appropriate protective equipment.

WARNING!

Damage to the respiratory tract and other organs due to escaping media.

- ▶ When sampling hazardous substances, follow the safety data sheet of the medium to be sampled.
- ▶ Wear the appropriate protective clothing.

NOTE!

Before using, the instrument has to be cleaned to the extent of your (validated) cleaning process.

- ▶ Measures of manual cleaning have to be carried out in the cleaning/desinfection area and under consideration of occupational health and safety.
 - ▶ Pay attention to the material resistance.
- ▶ Do not use any force when operating the instrument.

Layout and function

- ▶ The instrument is suitable for sampling small quantities of powders.
- ▶ The instrument is made of stainless steel and polyamide (PA) (push-button).

- ▶ Exchangeable volume heads allow sample quantities from 0.2 up to 10 ml (see accessories).
- ▶ The instrument is maintenance-free and suitable for steam sterilisation without the push-button (up to 121 °C).

Parts in contact with medium

- ▶ Stainless steel 1.4404/AISI 316 L

Point out any damage in the packaging to the freight forwarder/supplier immediately.

Assembly

CAUTION!

Openings with sharp edges on the instrument. Danger of cutting.

- ▶ Keep hands well away from openings with sharp edges.

CAUTION!

Moving parts on the instrument. Risk of crushing.

- ▶ Operate the push button carefully.

1. Keep hold of the tip ① and unscrew the push-button ② with the other hand.
2. Pull apart the instrument.
3. Screw in the desired volume tip (if necessary unscrew the mounted volume tip before).
4. Reassemble the instrument in reverse order.

Safe sampling

CAUTION!

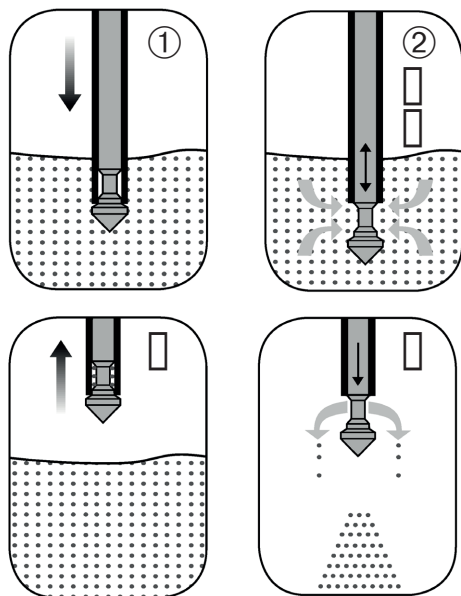
Pointed parts on the instrument. Risk of injury.

- ▶ Do not transport the instrument when the tip is not secured. Do not throw the instrument around.
- ▶ Do not pierce the instrument completely through the container.


NOTE!

The instrument is designed for manual operation.

- ▶ Do not hammer the instrument into the container by using a hammer.

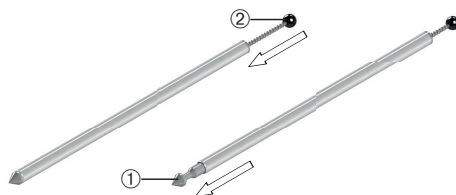1. Pierce the instrument into the medium to be sampled up to the desired depth.
2. At the desired depth, open the volume tip ① using the push-button ②.

3. The volume tip is filled with the sample.
4. Close the volume tip (release the push-button).
5. Pull the instrument out of the medium to be sampled.
6. Empty the sample into a suitable collecting container.

After taking the sample

1. Keep hold of the tip and unscrew the push-button with the other hand.
2. Pull apart the instrument.
3. Unscrew the tip.
4. Clean all parts.

Accessories

Type	Specification	Number	Item No.
MicroSampler heads	fits tube with 12 mm dia., 0,2 ml	1	5307-0102
MicroSampler heads	fits tube with 12 mm dia., 0,5 ml	1	5307-0105
MicroSampler heads	fits tube with 12 mm dia., 1,0 ml	1	5307-0110
MicroSampler heads	fits tube with 25 mm dia., 2,0 ml	1	5307-0202
MicroSampler heads	fits tube with 25 mm dia., 5,0 ml	1	5307-0205
MicroSampler heads	fits tube with 25 mm dia., 10,0 ml	1	5307-0210
Sample beaker with cover (PP, sterilisable)	125 ml	1	7103-0125

More accessories can be found at www.buerkle.de

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