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Directions for use | Instrucciones para el uso | Instruction d'emploi | Инструкция по пользованию

Art.-Nr. 5310-0150

Allschicht-Bohrer

All-layer drill

Perforador multicapa

Echantillonneur toutes couches

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General safety requirements

- ▶ Read these directions for use carefully before using.
- ▶ Keep these directions for use where they will be accessible for all users.
- ▶ Failure to comply with the directions for use shall render all warranty conditions null and void and may lead to personal injury and material damage.
- ▶ The instrument must only be used by qualified skilled employees.

Safety instructions before using

- ▶ Follow the manufacturer’s instructions and safety instructions of the medium you are sampling.
- ▶ Do not fill the instrument with potentially explosive substances.
- ▶ The instrument must only be used within the confines of applicable material resistance and for its designated purpose. Any other type of use will affect the warranty/guarantee.

Layout and function

The instrument is used for removing pourable powders and granulates. It is made of stainless steel. The instrument is maintenance-free and suitable for steam sterilisation (bis 121°C).

Parts in contact with medium

- ▶ Stainless steel 1.4404/AISI 316 L

Point out any damage in the packaging to the freight forwarder/supplier immediately.


Safe sampling

CAUTION!


Openings with sharp edges on the instrument.
Danger of cutting.

- ▶ Keep hands well away from openings with sharp edges.

1. Turn the upper handle ① on the instrument clockwise.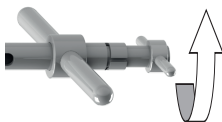


2. The instrument is closed and can pierce into the powder/granulate.
3. Screw the instrument with the aid of the lower handle ② to the desired depth.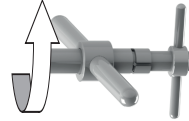


4. At the desired depth, turn the upper handle on the instrument counter clockwise.
5. The instrument opens and the sample is filled.
6. Turn the upper handle on the instrument clockwise.
7. The instrument is closed and can be removed from the powder/granulate.
8. Unscrew the instrument with the aid of the lower handle.
9. Turn the instrument around and empty the sample through the open inner tube into a suitable collecting container.

After taking the sample

1. Pull the inner tube all the way out by the upper handle ①.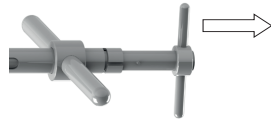

2. Unscrew the tip ②.

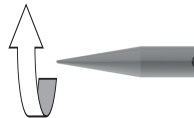3. Clean both tubes. Use the cleaning brushes (see accessories) for the inside of tubes.

Accessories

| Type | Specification | Number | Item No. |
|--|--------------------------------|--------|-----------|
| Cleaning brush | PVC/stainless steel, 100 cm | 1 | 5304-0100 |
| Control seal close-it | Printed, red, 95 x 95 mm | 1 | 5303-0012 |
| Control seal close-it | Printed, blue, 95 x 95 mm | 1 | 5303-0015 |
| Control seal close-it | Unprinted, white, 95 x 95 mm | 1 | 5303-0002 |
| Control seal close-it maxi | Printed, red, 150 x 150 mm | 1 | 5303-1012 |
| Control seal close-it maxi | Printed, blue, 150 x 150 mm | 1 | 5303-1015 |
| Control seal close-it maxi | Unprinted, white, 150 x 150 mm | 1 | 5303-1002 |
| Control seal close-it food & pharma | Printed, red, 95 x 95 mm | 1 | 5303-2017 |
| Control seal close-it food & pharma | Printed, blue, 95 x 95 mm | 1 | 5303-2021 |
| Control seal close-it food & pharma | Unprinted, white, 95 x 95 mm | 1 | 5303-2018 |
| Control seal close-it food & pharma maxi | Printed, red, 150 x 150 mm | 1 | 5303-3017 |
| Control seal close-it food & pharma maxi | Printed, blue, 150 x 150 mm | 1 | 5303-3021 |
| Control seal close-it food & pharma maxi | Unprinted, white, 150 x 150 mm | 1 | 5303-3018 |
| close-it tape | 95 mm x 50 m | 1 | 5303-0017 |
| close-it tape | 150 mm x 50 m | 1 | 5303-1017 |
| close-it tape food & pharma | 95 mm x 50 m | 1 | 5303-2019 |
| close-it tape food & pharma | 150 mm x 50 m | 1 | 5303-3019 |
| Tubes with screw cap, PP | 250 ml | 1 | 6214-0250 |
| Tubes with screw cap, PP | 600 ml | 1 | 6214-0600 |

More accessories can be found at www.buerkle.de

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