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Directions for use | Instrucciones para el uso | Instruction d'emploi | Инструкция по пользованию

Art.-Nr. 5315-0050 / 5315-0100 / 5315-0200

### Stechheber

Dip Sampler, plunging siphon

Pipeta

Siphon

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### General safety requirements

- ▶ Read these directions for use carefully before using.
- ▶ Keep these directions for use where they will be accessible for all users.
- ▶ Failure to comply with the directions for use shall render all warranty conditions null and void and may lead to personal injury and material damage.
- ▶ The instrument must only be used by qualified skilled employees.

### Safety instructions before using

- ▶ Follow the manufacturer's instructions and safety instructions of the medium you are sampling.
- ▶ When sampling hazardous substances, follow the technical rules on hazardous substances (in accordance with TRGS – (German) Technical Rules on Hazardous Substances) and wear appropriate protective equipment.
- ▶ The instrument must only be used within the confines of applicable material resistance and for its designated purpose. Any other type of use will affect the warranty/guarantee.
- ▶ Do not use any force when operating the instrument.

#### **WARNING!**

Damage to the respiratory tract and other organs due to escaping or outgasing media.

- ▶ When sampling hazardous substances, follow the safety data sheet of the medium to be sampled.
- ▶ Wear the appropriate protective clothing.

#### **NOTE!**

Before using, the instrument has to be cleaned to the extent of your (validated) cleaning process.

- ▶ Measures of manual cleaning have to be carried out in the cleaning/desinfection area and under consideration of occupational health and safety.
- ▶ Pay attention to the material resistance.

### Safety instructions during and after use

- ▶ The device can be used in the temperature range of 0 - 60 °C.
- ▶ Do not perform any lever movements during the application. The instrument may be deformed.

### Layout and function

- ▶ The instrument is made of PP.
- ▶ The instrument is suitable to take samples of liquids.
- ▶ The instrument is only suitable to take samples and not for continuous operation.
- ▶ The instrument is not suitable for the storage of samples.

### Parts in contact with medium

- ▶ PP
- ▶ HDPE

Point out any damage in the packaging to the freight forwarder/supplier immediately.

### Safe sampling

#### **WARNING!**

Bodily injury due to uncontrolled escaping liquid under pressure.

- ▶ Wear the appropriate protective clothing.

### Point sampling/soil sample

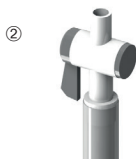
1. Close the tap ①.
2. Submerge the instrument into the medium to be sampled up to the desired depth / to the ground.
3. Open the tap ②. The instrument is filled with the liquid to be sampled.
4. Close the tap ①.
5. Remove the instrument from the medium to be sampled.
6. Turn the instrument over and empty the sample through the open tap into a suitable collecting container.

### All-layer sample

1. Open the tap ②.
2. Submerge the instrument into the medium to be sampled up to the desired depth / to the ground. Thereby, the instrument is filled with the liquid to be sampled.
3. Close the tap ①.
4. Remove the instrument from the medium to be sampled.
5. Turn the instrument over and empty the sample through the open tap into a suitable collecting container.

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