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Directions for use | Instrucciones para el uso | Instruction d'emploi | Инструкция по пользованию

Art.-Nr. 5326-0001 / 5326-0002

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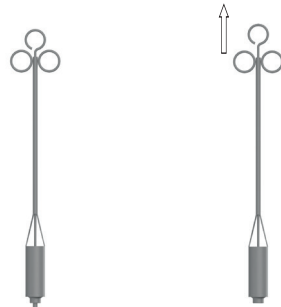
5. Gerät reinigen.

General safety requirements

- ▶ Read these directions for use carefully before using.
- ▶ Keep these directions for use where they will be accessible for all users.
- ▶ Failure to comply with the directions for use shall render all warranty conditions null and void and may lead to personal injury and material damage.
- ▶ The instrument must only be used by qualified skilled employees.

Safety instructions before using

- ▶ Follow the manufacturer's instructions and safety instructions of the medium you are sampling.
- ▶ When sampling hazardous substances, follow the technical rules on hazardous substances (in accordance with TRGS – (German) Technical Rules on Hazardous Substances) and wear appropriate protective equipment.
- ▶ The instrument must only be used within the confines of applicable material resistance and for its designated purpose. Any other type of use will affect the warranty/guarantee.
- ▶ Do not use any force when operating the instrument.

! WARNING!

Damage to the respiratory tract and other organs due to escaping or outgassing media.

- ▶ When sampling hazardous substances, follow the safety data sheet of the medium to be sampled.
- ▶ Wear the appropriate protective clothing.

NOTE!

Before using, the instrument has to be cleaned to the extent of your (validated) cleaning process.

- ▶ Measures of manual cleaning have to be carried out in the cleaning/desinfection area and under consideration of occupational health and safety.
- ▶ Pay attention to the material resistance.

Safety instructions during and after use

- ▶ The device can be used in the temperature range of 0 - 60 °C.
- ▶ Do not perform any lever movements during the application. The instrument may be deformed.
- ▶ The sampling area has to be free of disturbing factors and hazard sources (e.g. rotating elements).

Layout and function

- ▶ The instrument is made of stainless steel 1.4404 (AISI 316 L) / 1.4571 (AISI 316 Ti)
- ▶ The instrument is suitable to take samples of liquids and low-viscosity media without solid components.
- ▶ The instrument is only suitable to take samples and not for continuous operation.
- ▶ The instrument is not suitable for the storage of samples.
- ▶ The instrument is maintenance-free and suitable for steam sterilisation (up to 121°C).

Parts in contact with medium

- ▶ Stainless steel 1.4404 (AISI 316 L) / 1.4571 (AISI 316 Ti)

Point out any damage in the packaging to the freight forwarder/supplier immediately.

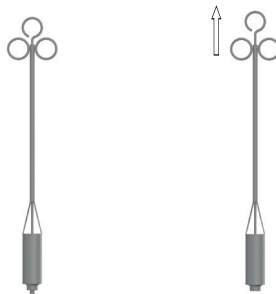
! WARNING!

Bodily injury due to uncontrolled escaping liquid.

- ▶ Check all parts for a firm fit and damages before use.
- ▶ Do not overturn the instrument when it is filled.
- ▶ Wear the appropriate protective clothing.

Safe sampling

1. Submerge the instrument into the medium to be sampled up to the desired depth.
2. The instrument is filled with the liquid to be sampled.
3. Remove the instrument from the medium to be sampled.
4. Open the valve ① using the thumb opening ② and empty the sample into a suitable sample container.

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