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Directions for use | Instrucciones para el uso | Instruction d'emploi | Инструкция по пользованию

Art.-Nr. 5340-1100

### Silo Probenehmer SiloPicker

Silo sampler SiloPicker

Muestreador Silo-picker

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### General safety requirements

- ▶ Read these directions for use carefully before using.
- ▶ Keep these directions for use where they will be accessible for all users.
- ▶ Failure to comply with the directions for use shall render all warranty conditions null and void and may lead to personal injury and material damage.
- ▶ The instrument must only be used by qualified skilled employees.

### Safety instructions before using

- ▶ Follow the manufacturer's instructions and safety instructions of the medium you are sampling.
- ▶ When sampling hazardous substances, follow the technical rules on hazardous substances (in accordance with TRGS – (German) Technical Rules on Hazardous Substances) and wear appropriate protective equipment.
- ▶ The instrument must only be used within the confines of applicable material resistance and for its designated purpose. Any other type of use will affect the warranty/guarantee.
- ▶ Do not use any force when operating the instrument.

#### **WARNING!**

Damage to the respiratory tract and other organs due to escaping or outgassing media.

- ▶ When sampling hazardous substances, follow the safety data sheet of the medium to be sampled.
- ▶ Wear the appropriate protective clothing.

#### **NOTE!**

Before using, the instrument has to be cleaned to the extent of your (validated) cleaning process.

- ▶ Measures of manual cleaning have to be carried out in the cleaning/desinfection area and under consideration of occupational health and safety.
- ▶ Pay attention to the material resistance.

### Safety instructions during and after use

- ▶ The device can be used in the temperature range of 0 - 60 °C.
- ▶ Pay attention to have sufficient space above during the application (minimum height: height of the sampler + 50 cm).

### Layout and function

- ▶ The instrument is made of stainless steel 1.4404/ AISI 316 L.
- ▶ The instrument is suitable to take bulk goods samples from silos.
- ▶ Using the extension rods (see accessories), samples can be taken from depths down to a maximum of 3.50 m.
- ▶ The insertion depth depends on the density of the bulk goods.
- ▶ The instrument is maintenance-free and suitable for steam sterilisation (up to 121°C).

### Parts in contact with medium

- ▶ Stainless steel 1.4404/AISI 316 L

### Scope of delivery

1 x	Sample container	①
1 x	Extension rod, 100 cm	②
1 x	Handle	③
1 x	Directions for use	

- ▶ Check delivery for completeness and transport damage.
- ▶ Point out any damage in the packaging to the freight forwarder/supplier immediately.

### Assembly

#### **CAUTION!**

- ▶ Moving parts on the instrument. Risk of crushing.
- ▶ Keep hands well away from moving parts.

1. Screw the handle ③ onto the extension rod ②.
2. If necessary, screw on additional extension rods (see accessories). Maximum permissible length: 3.50 m.
3. Screw the sample container ① onto the extension rod.

#### **NOTE!**

Possible loss of the sample container during use.

- ▶ Screw the sample container tightly onto the extension rod.
- ▶ Verify that the sample container is attached correctly before use.
- ▶ Do not rotate the instrument while plunging into the medium to be sampled.

**CAUTION!**

Bodily injury due to uncontrolled swinging around of the instrument and the extension rods.

- ▶ Pay attention to have sufficient space during the application.
- ▶ Disassemble the instrument for transport.

**Safe sampling****⚠ WARNING!**

Bodily injury due to accidental contact with electrical power. Danger of electric shock.

- ▶ Keep sufficient distance to any electrical power sources and supply lines.

**CAUTION!**

Pointed parts on the instrument. Risk of injury.

- ▶ Do not pierce the instrument completely through the container.
- ▶ Keep persons and body parts out of the danger area.

**CAUTION!**

Deformation of the extension rod. Slipping danger.

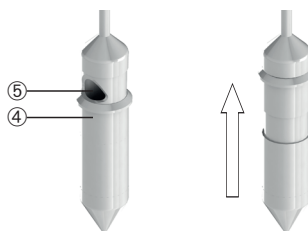
- ▶ Extend the rod only as long as necessary (maximum length 3.50 m).
- ▶ Do not plunge the instrument into highly compressed material.
- ▶ Do not perform any lever movements.

**CAUTION!**

Entanglement of the sampler by moving parts in the silo.

- ▶ Do not use the instrument, if moving parts in the silo are operating.

1. Pierce the instrument into the medium to be sampled up to the desired depth. Thereby, the sliding sleeve ④ closes the filling opening ⑤.

2. At the desired depth, pull back the instrument approx. 10 cm. Thereby, the sliding sleeve opens the filling opening.
3. Move the instrument gently up and down. Thereby, the sample container is filled with the medium to be sampled.
4. Pull the instrument out of the medium to be sampled.
5. Screw off the lower part of the sample container and empty the sample into a suitable collecting container.
6. Clean the instrument.

**Accessories**

Type	Specification	Number	Item No.
SiloPicker extension 50 cm	Stainless steel 1.4404/AISI 316 L	1	5380-0508
SiloPicker extension 100 cm	Stainless steel 1.4404/AISI 316 L	1	5380-1008
SiloPicker handle	Stainless steel 1.4404/AISI 316 L	1	0006-8112
SiloPicker sample container	Stainless steel 1.4404/AISI 316 L	1	5340-0100
Transport case	LxWxH 123x25x11 cm	1	3600-2120
Sample beaker with cover	PP, 125 ml	1	7103-0125
Sample container PFA	180 ml	1	6205-0180

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