**Bürkle GmbH**

Rheinauenstr. 5 | 79415 Bad Bellingen  
Telefon +49 (0) 76 35 6121-0  
Telefax +49 (0) 76 35 6121-31  
info@buerkle.de | www.buerkle.de

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### General safety requirements

- ▶ Read these directions for use carefully before using.
- ▶ Keep these directions for use where they will be accessible for all users.
- ▶ Failure to comply with the directions for use shall render all warranty conditions null and void and may lead to personal injury and material damage.
- ▶ The instrument must only be used by qualified skilled employees.

### Safety instructions before using

- ▶ Follow the manufacturer's instructions and safety instructions of the medium you are sampling.
- ▶ When sampling hazardous substances, follow the technical rules on hazardous substances (in accordance with TRGS – (German) Technical Rules on Hazardous Substances) and wear appropriate protective equipment.
- ▶ The instrument must only be used within the confines of applicable material resistance and for its designated purpose. Any other type of use will affect the warranty/guarantee.
- ▶ Do not use any force when operating the instrument

#### NOTE!

Before using, the instrument has to be cleaned to the extent of your (validated) cleaning process.

- ▶ Measures of manual cleaning have to be carried out in the cleaning/desinfection area and under consideration of occupational health and safety.
- ▶ Pay attention to the material resistance.

### Safety instructions during and after use

- ▶ The instrument is not suitable for the use with wood, stone and metal.
- ▶ Do not use the instrument on live humans or animals.
- ▶ Wear the appropriate protective clothing.

### Layout and function

- ▶ The instrument is made of stainless steel 1.4034 (AISI 420).
- ▶ The instrument is suitable to take samples of semi-solid materials in the food and beverage industry using hand-held drills or cordless screwdrivers.
- ▶ The instrument is only suitable to take samples and not for continuous operation.
- ▶ The instrument is maintenance-free and suitable for

steam sterilisation (up to 121°C) for sterilization by flame treatment.

- ▶ If necessary, the cutting edge can be resharpened.

### Parts in contact with medium

- ▶ Stainless steel 1.4034 (AISI 420).

Point out any damage in the packaging to the freight forwarder/supplier immediately.

### Assembly

1. Mount the drill head  on a hand-held drill or cordless screwdriver.

#### NOTE!

Verify that the borer is attached correctly before use.

### Safe sampling

#### WARNING!

Bodily injury due to accidental contact with electrical power. Danger of electric shock.

- ▶ Keep the cable of the drill out of the cutting area.

#### WARNING!

Sharp parts on the instrument. Risk of injury.

- ▶ Keep hands well away from the cutting area.

#### WARNING!

Bodily injury due to drilling through the sampling material without adequate securing.

- ▶ Use a suitable base (e.g. cutting board made of plastic or wood).
- ▶ Use a suitable locking mechanism or secure the sampling material against twisting and shifting.
- ▶ Only place the instrument vertically onto the sampling material.

#### WARNING!

Bodily injury due to entanglement of body parts, hairs or clothes.

- ▶ Keep hands well away from the cutting area.
- ▶ Use a hairnet if necessary.
- ▶ Ensure to wear tight-fitting clothes.

1. Place the drill head ① onto the sampling material.
2. Start boring slowly until a slight trace is incised in the sampling material.
3. Pull the drill head ① out of the sampling material.
4. Remove the sample carefully from the sampling material.
5. Clean the instrument.

**CAUTION!**

Sharp parts on the instrument. Risk of injury.

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