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Directions for use | Instrucciones para el uso | Instruction d'emploi | Инструкция по пользованию

Art.-Nr. 5389-0010 / 5389-0100 / 5389-0250 / 5389-0500 / 5389-1000

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PTFE Scoop

Cuchara de PTFE

Puisard PTFE

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### General safety requirements

- ▶ Read these directions for use carefully before using.
- ▶ Keep these directions for use where they will be accessible for all users.
- ▶ Failure to comply with the directions for use shall render all warranty conditions null and void and may lead to personal injury and material damage.
- ▶ The instrument must only be used by qualified skilled employees.

### Safety instructions before using

- ▶ Follow the manufacturer's instructions and safety instructions of the medium you are sampling.
- ▶ When sampling hazardous substances, follow the technical rules on hazardous substances (in accordance with TRGS – (German) Technical Rules on Hazardous Substances) and wear appropriate protective equipment.
- ▶ The instrument must only be used within the confines of applicable material resistance and for its designated purpose. Any other type of use will affect the warranty/guarantee.

#### **WARNING!**

Damage to the respiratory tract and other organs due to escaping or outgassing media.

- ▶ When sampling hazardous substances, follow the safety data sheet of the medium to be sampled.
- ▶ Wear the appropriate protective clothing.
- ▶ The instrument is only suitable to take samples and not for continuous operation.
- ▶ Only use the instrument under adequate weather conditions.
- ▶ Do not use the instrument in frozen watercourses.
- ▶ Operate the instrument only with the human power of a person.
- ▶ Secure the sampling area before sampling (pay attention to humans, objects, ships, moving parts etc.).
- ▶ Do not use any force when operating the instrument.

#### **NOTE!**

Before using, the instrument has to be cleaned to the extent of your (validated) cleaning process.

- ▶ Measures of manual cleaning have to be carried out in the cleaning/desinfection area and under consideration of occupational health and safety.
- ▶ Pay attention to the material resistance.

### Safety instructions during and after use

- ▶ The instrument is temperature-resistant from -270 to +270 °C.
- ▶ Pay attention to have sufficient space during use: Radius of the length of the scoop + 50 cm.
- ▶ Only perform lifting, lowering and slight lateral movements during use.

### Layout and function

- ▶ The instrument is made of PTFE.
- ▶ The instrument is suitable to take samples of liquids and low-viscous substances.
- ▶ The instrument is not suitable to take samples of bulk goods.
- ▶ The 60 cm long rod can be extended to 120 cm by adding another component (see accessories).
- ▶ The scoop PTFE with 10 ml has a strong rod with a length of 155 mm.
- ▶ The instrument is maintenance-free.

### Parts in contact with medium

- ▶ PTFE

### Assembly

If necessary mount the extension rod (see accessories).

### Safe sampling

#### **WARNING!**

Risk of slipping and falling due to wet or smooth floor.

- ▶ Ensure to have enough support and stability.

#### **WARNING!**

Breakage during sampling.

- ▶ Do not overload the scoop.
- ▶ Perform a visual check for material fatigue (cracks etc.) before sampling.
- ▶ Ensure that the scoop is attached correctly before use.

#### **CAUTION!**

Deformation or breakage of the rod.

- ▶ Do not perform any lever movements.

**CAUTION!**

Bodily injury due to uncontrolled swinging around of the instrument.

- ▶ Pay attention to have sufficient space during use: Radius of the length of the scoop + 50 cm.

1. Take the sample using the scoop.
2. Empty the sample into a suitable collecting container.
3. Clean the instrument.

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