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Directions for use | Instrucciones para el uso | Instruction d'emploi | Инструкция по пользованию

Art.-Nr. 5601-0130

Fassverschraubung gasdicht

Gastight barrel connector

Abrazadera de rosca estanca al gas

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General safety requirements

- ▶ Read these directions for use carefully before using.
- ▶ Keep these directions for use where they will be accessible for all users.
- ▶ Failure to comply with the directions for use shall render all warranty conditions null and void and may lead to personal injury and material damage.
- ▶ The instrument must only be used by qualified skilled employees.
- ▶ The instrument is designed for use in industry, laboratories and science. The instrument is not suitable for any application in private or public spaces.

Safety instructions before using

- ▶ Follow the manufacturer's instructions and safety instructions of the medium you are filling.
- ▶ Damage to the respiratory tract and other organs due to escaping or off-gassing media. Fill only in ventilated rooms or under a suitable exhaust hood.

WARNING!

Possible sparking due to electrostatic discharge. Danger of explosion if combustible liquids are used.

- ▶ Thread sealing tape influences the electrical conductivity. The barrel connector is equipped with a sealing ring. No sealing tape necessary.
 - ▶ Connect the earthing terminal before use (see accessories).
 - ▶ Wear the necessary protective clothing.
- ▶ The instrument must only be used within the confines of applicable material resistance and for its designated purpose. Any other type of use will affect the warranty/guarantee.

Safety instructions during and after use

- ▶ The instrument is suitable at a temperature range from 0 to 60 °C.
- ▶ The instrument is not suitable for cleaning in a dishwasher.

Layout and function

- ▶ The instrument is made of Nickel-plated brass, the seals made of FKM.
- ▶ The instrument is to attach manual pumps (immersion tube Ø 32 mm) in vertical position in barrels with connection thread R 2".

- ▶ The instrument is suitable for the gas-tight filling of liquids and slightly viscous media in conjunction with a manual pump (see accessories). A one-sided air aeration is given by a non return-valve.

WARNING!

Possible escaping of liquid and gas due to excess pressure in the barrel. Bodily injury due to escaping gas and liquid.

- ▶ Avoid excessive excess pressure. (Do not expose the barrel to direct solar irradiation or variations in temperature. Pay attention to chemical resistance.)
- ▶ Ensure adequate air ventilation.
- ▶ Wear the necessary protective clothing.

NOTE!

Use a suitable pump to ensure gas-tight filling.

- ▶ It is recommended to use a pump that can be hermetically sealed by shut-off valves (see accessories).
- ▶ A gas-shuttle pipe can be connected to non return valve ⑤ by a tubing connector (thread connection 1/8").

NOTE!

Non return valve ⑤ is set to the correct partial vacuum in the factory. Do not adjust it.

Scope of delivery

1 x	Barrel connector	①
1 x	Allen wrench	②

- ▶ Check delivery for completeness and transport damage.
- ▶ Point out any damage in the packaging to the freight forwarder/supplier immediately.

Assembly

1. Screw on the barrel connector ① hand-tight and air-tight (R 2" steel barrel thread).

NOTE!

Regularly check seal for connection and leak tightness.

NOTE!

Check correct connection.

- ▶ Screw on barrel connector ① only hand tight and without force.

2. Slightly moisten immersion tube and insert it in the barrel connector ①.

NOTE!

Ensure to use a fitting immersion tube (Immersion tube Ø 32 mm) (see accessories).

3. Slide the pump until it has reached the required immersion depth.

4. Remove allen wrench ② from magnet holder ③ and tighten up the fixing screw ④.
5. Put the allen wrench ② back to magnet holder ③.

Dismounting

1. Remove allen wrench ② from magnet holder ③ and loosen the fixing screw ④.
2. Put the allen wrench ② back to magnet holder ③.
3. Carefully remove the pump from the barrel connector ①.
4. Unscrew the barrel connector ①.
5. Rinse the barrel connector ① with a pH-neutral fluid.

Accessories

Type	Specification	Number	Item No.
Stainless steel barrel pump	Discharge tube, immersion depth 36 cm	1	5601-0400
Stainless steel barrel pump	Discharge tube, immersion depth 57 cm	1	5601-0500
Stainless steel barrel pump	Discharge tube, immersion depth 91 cm	1	5601-0800
Stainless steel barrel pump	Discharge hose & Stopcock, immersion depth 36 cm	1	5601-0401
Stainless steel barrel pump	Discharge hose & Stopcock, immersion depth 57 cm	1	5601-0501
Stainless steel barrel pump	Discharge hose & Stopcock, immersion depth 91 cm	1	5601-0801
Anti-static set	To prevent electrostatic charging	1	5602-1000

More accessories can be found at www.buerkle.de

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