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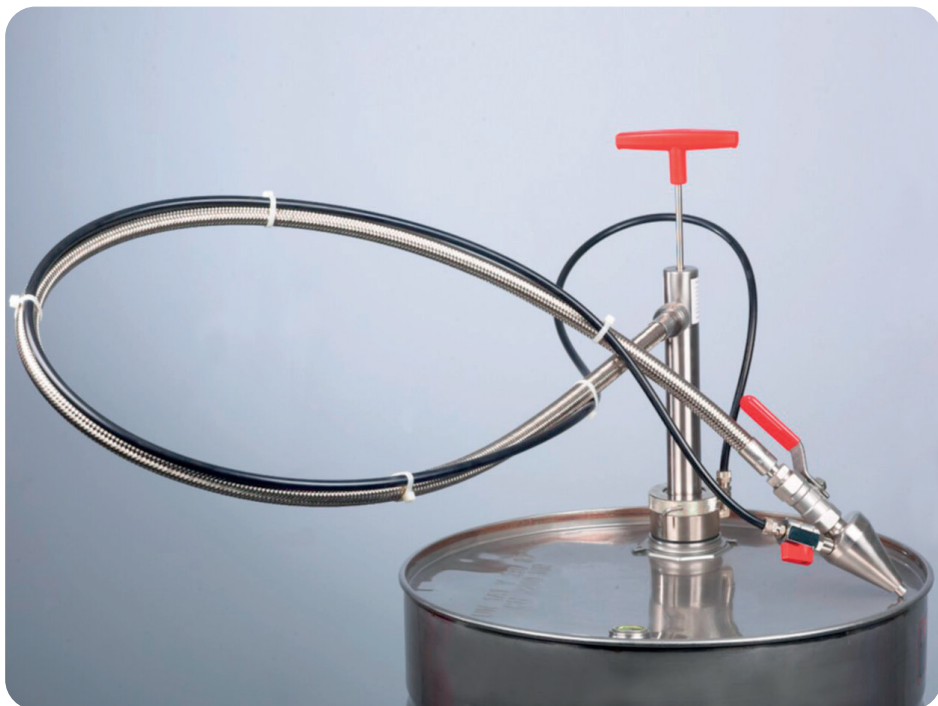
Directions for use | Instrucciones para el uso | Instruction d'emploi | Инструкция по пользованию

Art.-Nr. 5611-1090 / 5611-0090

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## General safety requirements

- ▶ Read these directions for use carefully before using.
- ▶ Keep these directions for use where they will be accessible for all users.
- ▶ Failure to comply with the directions for use shall render all warranty conditions null and void and may lead to personal injury and material damage.
- ▶ The instrument must only be used by qualified skilled employees.
- ▶ The instrument is designed for use in industry, laboratories and science. The instrument is not suitable for any application in private or public spaces.

## Safety instructions before using

- ▶ Follow the manufacturer's instructions and safety instructions of the medium you are filling.
- ▶ Damage to the respiratory tract and other organs due to escaping or off-gassing media. Fill only in ventilated rooms or under a suitable exhaust hood. Wear appropriate protective clothing.
- ▶ When working with explosive substances, follow the technical rules for combustible liquids (in accordance with TRBF – (German) Technical Rules on Combustible Liquids) and wear appropriate protective equipment.

### **WARNING!**

Possible sparking due to electrostatic discharge. Danger of explosion if combustible liquids are used.

- ▶ Connect the earthing terminal before use (see accessories).
- ▶ Wear the necessary protective clothing.
- ▶ The instrument must only be used within the confines of applicable material resistance and for its designated purpose. Any other type of use will affect the warranty/guarantee.
- ▶ Not suitable for acetone.
- ▶ Do not use any force when operating the instrument.

### **CAUTION!**

Escaping liquids due to falling over or tipping of the instrument. Bodily injury due to escaping liquid and damage to the instrument.

- ▶ Ensure sufficient stability of the instrument.
- ▶ Shut down and secure the instrument immediately.
- ▶ Contact the dealer or manufacturer.

## Safety instructions during and after use

- ▶ The instrument is suitable for indoor use and manual operation.
- ▶ Pump with manual operation. Not suitable for continuous operation.
- ▶ Pay attention to have sufficient space above during the filling process (minimum height: height of the maximum extended piston + 20 cm).
- ▶ The instrument is not suitable for cleaning in a dishwasher.

### **NOTE!**

Pump and valves can be damaged or plugged by dirt particles.

- ▶ Do not fill off any polluted media.
- ▶ Rinse the pump with a pH-neutral medium.

## Layout and function

- ▶ The instrument is made of conductive material (stainless steel (1.4301), nickel plated brass, PA electrically conductive) and PTFE. It is suitable to fill off combustible liquids.
- ▶ The instrument is suitable for filling from containers with a capacity of 60 - 1000 l into smaller containers.
- ▶ The instrument is maintenance-free and suitable at a temperature range from 0 to 40 °C.

Point out any damage in the packaging to the freight forwarder/supplier immediately.

### **WARNING!**

Possible electrostatic discharge. Danger of explosion if combustible liquids are used.

- ▶ Connect the earthing terminal before use (see accessories).
- ▶ Wear the necessary protective clothing.

## Assembly

### **WARNING!**

Some liquid will continue to run. Bodily injury due to escaping liquid.

- ▶ Close the shut-off valves ③ on the discharge hose ① and the gas-shuttle pipe ② before mounting.

1. Guide the barrel screw connection ⑤ and riser pipe ④ into the container and screw it together finger-tight.
2. If necessary adjust the riser pipe.

**NOTE!**

- Ensure to use a fitting thread (see accessories).  
 ▶ Screw the thread by hand and without any force.

3. Screw the discharge hose onto the connecting thread of the pump ⑥.
4. Attach the gas-shuttle pipe to the thread of the barrel screw joint ⑦ and screw it together finger-tight using the cap nut.
5. Attach the gas-shuttle pipe to the thread of the discharge hose ⑧ and screw it together finger-tight using the cap nut.

**NOTE!**

- ▶ Check the hose connections and seals regularly for position and tightness.

**Safe filling****CAUTION!**

Liquid entering the instrument. Bodily injury due to escaping liquid and damage to the instrument.

- ▶ Do not overfill the filling container.  
 ▶ Do not fill off when the shut-off valves are closed.
1. Open the shut-off valves on the discharge hose and the gas-shuttle pipe.
  2. Press the filling cone ⑨ into the opening of the filling container to form an airtight seal. The filling cone has to be pressed onto the container during the whole filling process.

**NOTE!**

- Only filling containers with round shaped openings and a diameter of 18 to 49 mm can be used.

3. Activate the piston rod ⑩ by using the handle and fill off the liquid. Thereby, the displaced air is returned from the filling container to the original container via the gas-shuttle pipe.

**Completing filling and dismantling**

1. After the last pump stroke, leave the piston rod in the lower position and insert it into the pump up to the stop.
2. Close both shut-off valves.
3. To prevent any liquid continue to run (siphoning principle), fix the discharge hose above the liquid level in the original container.

**! WARNING!**

Some liquid will continue to run. Bodily injury due to escaping liquid.

- ▶ Close the shut-off valves on the discharge tube and the gas-shuttle pipe before dismantling.  
 ▶ Raise the discharge hose and allow the liquid to run back into the container.
4. Carefully remove the pump from the original container and allow the remaining liquid to drain into the filling container.
  5. Allow the remaining vapours to escape out of the instrument and rinse the instrument with a pH-neutral liquid.

**Accessories**

Type	Specification	Number	Item No.
Antistatic set	To prevent electrostatic charging	1	5602-1000
SafetyBox	Incl. safety goggles and protective gloves	1	2507-9000

More accessories can be found at [www.buerkle.de](http://www.buerkle.de)

**Troubleshooting**

Problem	To eliminate the error
Liquid is not pumped.	Bottom valve is polluted. Clean the pump.
	The telescopic riser pipe does not extend into the liquid being filled. Open the shut-off valves.
	If there is only little residual liquid in the container, the pumping must be carried out for a longer time.

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