Bürkle GmbH

Rheinauenstr. 5 | 79415 Bad Bellingen
Telefon +49 7635 6121-0
info@buerkle.de
www.buerkle.de

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General safety requirements

- ▶ Read these directions for use carefully before using.
- ▶ Keep these directions for use where they will be accessible for all users.
- ▶ Failure to comply with the directions for use shall render all warranty conditions null and void and may lead to personal injury and material damage.
- ▶ The instrument must only be used by qualified skilled employees.

Safety instructions

- ▶ Follow the manufacturer's instructions and safety instructions of the medium you are filling.
- ▶ Damage to the respiratory tract and other organs due to off-gassing media. Fill only in well ventilated rooms or under a suitable exhaust hood.
- ▶ The instrument must only be used within the confines of applicable material resistance and for its designated purpose. Any other type of use will affect the warranty/guarantee.

⚠ WARNING!

Possible sparking due to electrostatic discharge. Danger of explosion if combustible liquids are used.

- ▶ Connect the earthing cable to the Connection (M6 thread) ③ before use.
- ▶ Wear the necessary protective clothing.

CAUTION!

Possible overturning of the drum due to insufficient securing.

- ▶ Make certain the drum is positioned with adequate stability.
- ▶ Take appropriate measures to prevent the drum from tipping over.

Layout and function

- ▶ The instrument is made of stainless steel V2A.
- ▶ The drum holder is suitable for placing filling containers (load capacity: 10 kg).
- ▶ Can be used for plastic or stainless steel drums with at least 600 mm, 10 mm rim height and maximum capacity of 60 litres to 220 litres.
- ▶ The storage area with adjustable height is 240 x 240 mm.
- ▶ When used with discharge hose the maximum height of the target container is 470 mm.

NOTE!

The holder can be used for Bürkle pumps with discharge tube or discharge hose (see accessories).

- ▶ The instrument is suitable at a temperature range from 0 to 60 °C.
- ▶ Constructional changes of the instrument or the accessories aren't allowed.

Point out any damage in the packaging to the freight forwarder/supplier immediately.

Assembly and safe usage

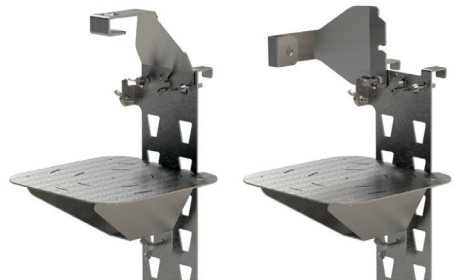
1. Hang the drum holder ① on the drum edge with the hook ②.

CAUTION!

Squeezing, cutting the fingers and body parts possible when mounting the holder.

- ▶ Mount the holder carefully.

2. Fix the drum holder ① in place with clamping screws ③.
3. Attach the storage area ④ at the desired height and by means of adjusting screws ⑤ if necessary put it in a horizontal position.
4. When used with discharge tube, fold away the Fastening for discharge hose ⑥. When used with discharge hose, straighten the Fastening for discharge hose ⑥ and, if necessary, fix it with the wing screw ⑦ (see figure). Insert the discharge hose.

With discharge hose With discharge tube

5. If necessary connect the earthing cable to the connection for earthing cable (M6 thread) ③.
6. Place the filling container under the discharge tube/hose.

Accessories

Type	Specification	Number	Item No.
PTFE barrel pump Ultrapure - discharge tube	Immersion depth 60 cm	1	5606-0600
PTFE barrel pump Ultrapure - discharge tube	Immersion depth 95 cm	1	5606-1000
PP barrel pump - discharge tube	Immersion depth 65 cm	1	5600-0651
PP barrel pump - discharge tube	Immersion depth 80 cm	1	5600-0801
PP barrel pump - discharge tube	Immersion depth 100 cm	1	5600-1001
PP barrel pump - discharge tube	Immersion depth 125 cm	1	5600-1251
PP barrel pump - discharge tube	Immersion depth 150 cm	1	5600-1501
Stainless steel barrel pump - discharge tube	Immersion depth 91 cm	1	5601-0800
Stainless steel barrel pump - discharge hose/stopcock	Immersion depth 91 cm	1	5601-0801
Solvent pump foot operated - discharge tube	Immersion depth 95 cm	1	5603-1100
Solvent pump foot operated - discharge hose/stopcock	Immersion depth 95 cm	1	5603-3000
Withdrawal system for solvents with discharge tube	Immersion depth 95 cm	1	5603-5003
Withdrawal system for solvents with discharge hose	Immersion depth 95 cm	1	5603-5004

More accessories can be found at www.buerkle.de

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