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Directions for use | Instrucciones para el uso | Instruction d'emploi | Инструкция по пользованию

Art.-Nr. 7118-2000

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General safety requirements

- ▶ Read these directions for use carefully before using.
- ▶ Keep these directions for use where they will be accessible for all users.
- ▶ Failure to comply with the directions for use shall render all warranty conditions null and void and may lead to personal injury and material damage.
- ▶ The instrument must only be used by qualified skilled employees.

Safety instructions before using

- ▶ Follow the manufacturer's instructions and safety instructions of the medium.
- ▶ Damage to the respiratory tract and other organs due to off-gassing media. Use only in ventilated rooms or under a suitable exhaust hood.
- ▶ The instrument must only be used within the confines of applicable material resistance and for its designated purpose. Any other type of use will affect the warranty/guarantee.
- ▶ The instrument is suitable for indoor use.
- ▶ The instrument is designed for stationary use and is not suitable for the transport of funnels.

Safety instructions during use

WARNING!

Danger due to unsuitable working area. Danger of tipping over.

- ▶ Bright, dry and clean working place at table height without obstructions.
 - ▶ Horizontal stand of the frame with sufficient space to fill the funnels.
- ▶ The instrument has to be mounted on a horizontal, even surface.
 - ▶ The device can be used in the temperature range of 0 - 40 °C.

Layout and function

- ▶ The instrument consists of a foot and two holders for sedimentation funnels made of polypropylen (PP) and a support rod made of chrome-plated steel.
- ▶ The instrument is suitable to hold sedimentation funnels during the sedimentation.
- ▶ The instrument is suitable for funnels with a diameter of at least 90 mm and a maximum capacity of 1 litre.
- ▶ The instrument is maintenance-free.


Scope of delivery

1 x	Support rod, Ø 16 mm, chrome-plated steel	①
1 x	Foot with two deposit spaces, PP	②
2 x	Holder for sedimentation funnels with magnets (each consisting of two funnel holders with different diameters), PP	③ ④
1 x	Directions for use	


- ▶ Check delivery for completeness and transport damage.
- ▶ Point out any damage in the packaging to the freight forwarder/supplier immediately.

Assembly

1. Screw off the screw from the support rod ①.
2. Insert the support rod from above into the foot ②.
3. Fasten the support rod with the aid of the screw from below (by rotating the support rod).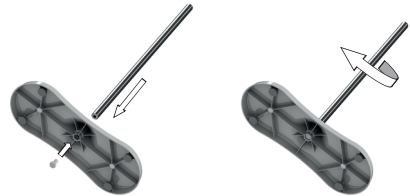

4. Plug the funnel holders ③ ④ onto the support rod and move them to the desired height.
5. Position the holders above the deposit spaces ⑤.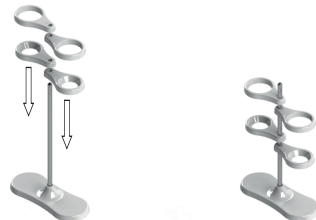


Safe usage**⚠ WARNING!**

Bodily injury due to tipping over of the instrument.

- ▶ Position the funnels only above the deposit spaces (see diagram).
- ▶ Maximum load per funnel holder: 2 kg.

1. Position the funnel holders above the deposit spaces.
2. Insert the funnels into the funnel holders.
3. Fill in the medium and run the sedimentation.

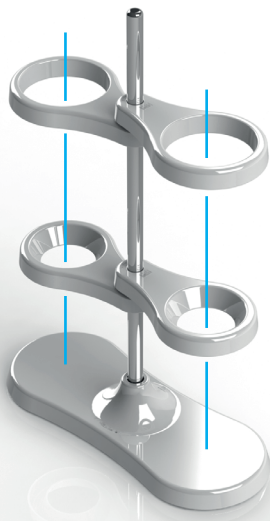Correct position of the funnel holders

Wrong position of the funnel holders -Danger of tipping over!

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