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Directions for use | Instrucciones para el uso | Instruction d'emploi | Инструкция по пользованию

Art.-Nr. 7118-2001

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### General safety requirements

- ▶ Read these directions for use carefully before using.
- ▶ Keep these directions for use where they will be accessible for all users.
- ▶ Failure to comply with the directions for use shall render all warranty conditions null and void and may lead to personal injury and material damage.
- ▶ The instrument must only be used by qualified skilled employees.

### Safety instructions before using

- ▶ Follow the manufacturer's instructions and safety instructions of the medium.
- ▶ Damage to the respiratory tract and other organs due to off-gassing media. Use only in ventilated rooms or under a suitable exhaust hood.
- ▶ The instrument must only be used within the confines of applicable material resistance and for its designated purpose. Any other type of use will affect the warranty/guarantee.
- ▶ The instrument is suitable for indoor use.
- ▶ The instrument is designed for stationary use and is not suitable for the transport of funnels.

### Safety instructions during use

#### **WARNING!**

Danger due to unsuitable working area. Danger of tipping over.

- ▶ Bright, dry and clean working place at table height without obstructions.
  - ▶ Horizontal stand of the frame with sufficient space to fill the funnels.
- ▶ The instrument has to be mounted on a horizontal, even surface.
  - ▶ The device can be used in the temperature range of 0 - 40 °C.

### Layout and function

- ▶ The instrument is suitable to hold sedimentation funnels during the sedimentation.
- ▶ The instrument is suitable for funnels with a diameter of at least 90 mm and a maximum capacity of 1 litre.
- ▶ The instrument is maintenance-free.


### Scope of delivery

1 x	Holder for sedimentation funnels with magnets (each consisting of two funnel holders with different diameters), PP	③ ④
1 x	Directions for use	

- ▶ Check delivery for completeness and transport damage.
- ▶ Point out any damage in the packaging to the freight forwarder/supplier immediately.

### Assembly

1. Plug the funnel holders ③ ④ onto the support rod and move them to the desired height.
2. Position the holders above the deposit spaces ⑤.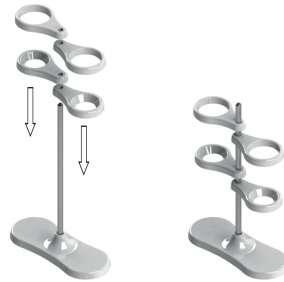


### Safe usage


#### **WARNING!**

Bodily injury due to tipping over of the instrument.

- ▶ Position the funnels only above the deposit spaces (see diagram).
  - ▶ Maximum load per funnel holder: 2 kg.
1. Position the funnel holders above the deposit spaces.
  2. Insert the funnels into the funnel holders.
  3. Fill in the medium and run the sedimentation.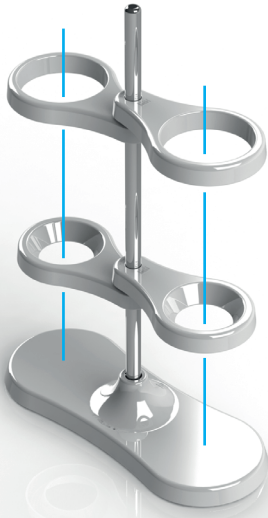

**Correct position of the funnel holders**

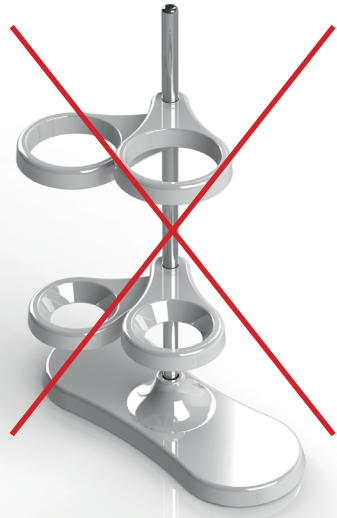**Wrong position of the funnel holders -Danger of tipping over!**

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