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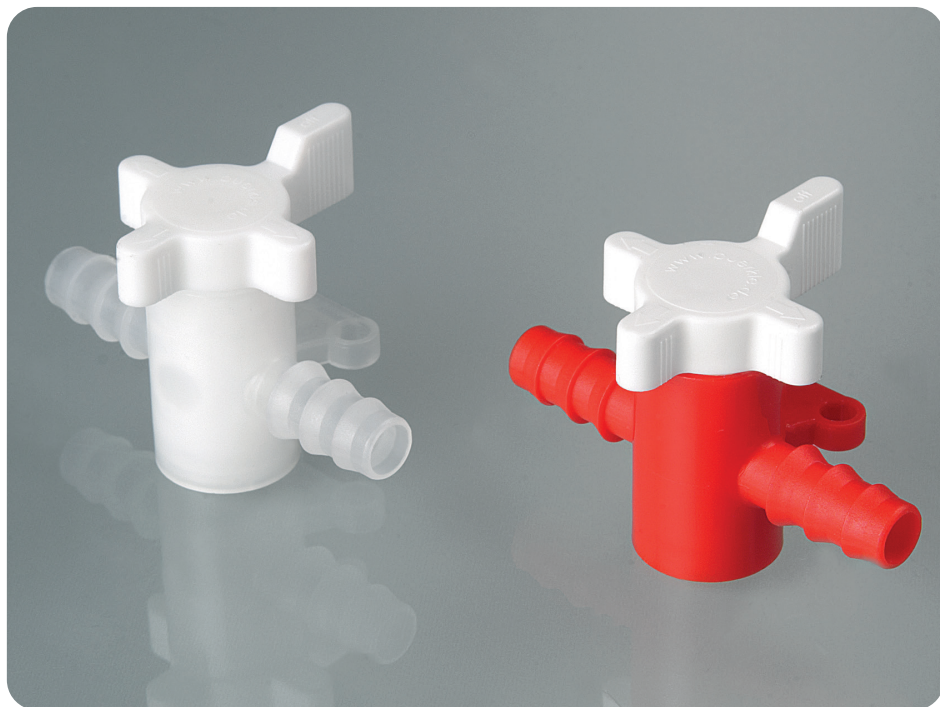
Directions for use | Instrucciones para el uso | Instruction d'emploi | Инструкция по пользованию

Art.-Nr. 8605-0060 / 8605-0080 / 8605-0100 / 8605-0120 / 8607-0060 /
8607-0080 / 8607-0100 / 8607-0120

Zweiwegeventil

Two-way valves

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General safety requirements

- ▶ Read these directions for use carefully before using.
- ▶ Keep these directions for use where they will be accessible for all users.
- ▶ Failure to comply with the directions for use shall render all warranty conditions null and void and may lead to personal injury and material damage.
- ▶ The instrument must only be used by qualified skilled employees.

Safety instructions

- ▶ Follow the manufacturer's instructions and safety instructions of the medium to be transferred.
- ▶ The instrument must only be used within the confines of applicable material resistance and for its designated purpose. Any other type of use will affect the warranty/guarantee.
- ▶ The instrument is suitable for indoor use and manual operation. Do not expose the device to any direct sunlight.
- ▶ The instrument is not suitable for cleaning in a dishwasher.

NOTE!

Before using, the instrument has to be cleaned to the extent of your (validated) cleaning process.

- ▶ Measures of manual cleaning have to be carried out in the cleaning/desinfection area and under consideration of occupational health and safety.
- ▶ Pay attention to the material resistance.

Layout and function

- ▶ The device is made of PP/PE and PVDF.
- ▶ The device is suitable for releasing or blocking the flow of gases and liquids.

- ▶ The device is suitable for liquids (pay attention to chemical resistance) and hazard-free gases (air).
- ▶ The device can be used in the temperature range of 5 to 40 °C.
- ▶ The device is pressure-resistant up to max. 1 bar.

Point out any damage in the packaging to the freight forwarder/supplier immediately.

Assembly and safe usage

1. If necessary, fasten the two-way valve ① using the mounting plate ②.

NOTE!

The installation of the device is only allowed by trained specialists.

- ▶ The user is responsible for the correct mounting in connection with the substrate condition.
- ▶ Drill hole diameter is 4.2 mm (fixing material not included).


2. Attach the hoses to the threads ③.

NOTE!

If necessary, fasten hoses with hose clips (see accessories).

- ▶ Do not overtighten hose clips.

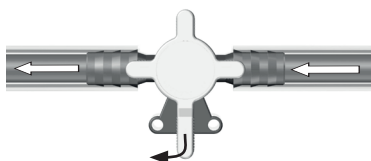3. Turn the plug ④ to release the flow.

CAUTION!

Possible danger of breakage during operation.

- ▶ Apply only normal manual force.

CAUTION!

Spaces on the device. Risk of crushing.

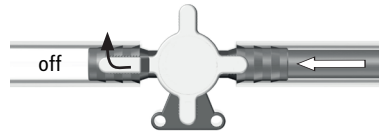
- ▶ Be careful for spaces when activating the plug.

! WARNING!

Danger of bodily injury due to liquid under high pressure escaping.

- ▶ The pressure must not exceed 1 bar.
- ▶ Wear the necessary protective clothing.

4. Turn the plug ④ further to block the flow.

! WARNING!

Bodily injury due to escaping liquid.

- ▶ Do not remove the device from filled hoses.
- ▶ Empty hoses and unit completely before disassembly.

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