Bürkle GmbH

Rheinauenstr. 5 | 79415 Bad Bellingen
Telefon +49 7635 6121-0
info@buerkle.de
www.buerkle.de

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General safety requirements

- ▶ Read these directions for use carefully before using.
- ▶ Keep these directions for use where they will be accessible for all users.
- ▶ Failure to comply with the directions for use shall render all warranty conditions null and void and may lead to personal injury and material damage.
- ▶ The instrument must only be used by qualified skilled employees.

Safety instructions before using

- ▶ Follow the manufacturer's instructions and safety instructions of the medium you are filling.
- ▶ Damage to the respiratory tract and other organs due to off-gassing media. Use only in ventilated rooms or under a suitable exhaust hood.
- ▶ The instrument must only be used within the confines of applicable material resistance and for its designated purpose. Any other type of use will affect the warranty/guarantee.
- ▶ The instrument is suitable for indoor use.
- ▶ The instrument is designed for stationary use and is not suitable for the transport of funnels.

Safety instructions during use

WARNING!

Danger due to unsuitable working area. Danger of tipping over.

- ▶ Bright, dry and clean working place at table height without obstructions.
 - ▶ Horizontal stand of the frame with sufficient space to fill the funnels.
- ▶ The instrument has to be mounted on a horizontal, even surface.
 - ▶ The device can be used in the temperature range of 0 - 40 °C.
 - ▶ The instrument may only be used for funnels.

Layout and function

- ▶ The instrument consists of a foot and two funnel holders made of polypropylen (PP), two reducing inserts made of PVC and a support rod made of chrome-plated steel.
- ▶ The instrument is suitable to hold funnels. It simplifies filling and filtering.

- ▶ The instrument is suitable for funnels with an upper outer diameter from 40 to a maximum of 180 mm (use the reducing inserts for funnels smaller than 80 mm) and a maximum capacity of 1 litre.
- ▶ The instrument is maintenance-free.


Scope of delivery

1 x	Support rod, Ø 16 mm, chrome-plated steel	①
1 x	Foot with two deposit spaces, PP	②
2 x	Funnel holders with magnets, PP	③
2 x	Reducing inserts for small funnels, PVC	④
1 x	Directions for use	


- ▶ Check delivery for completeness and transport damage.
- ▶ Point out any damage in the packaging to the freight forwarder/supplier immediately.

Assembly

1. Screw off the screw from the support rod ①.
2. Insert the support rod from above into the foot ②.
3. Fasten the support rod with the aid of the screw from below (by rotating the support rod).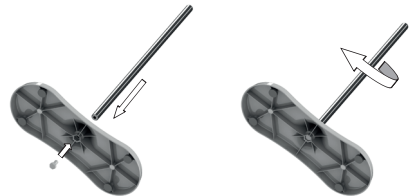

4. Plug the funnel holders ③ onto the support rod and move them to the desired height. Position the holders above the deposit spaces ⑤.

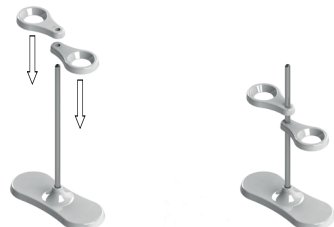5. If necessary, place the reducing inserts ④ into the funnel holders (when using small funnels).


Safe usage

WARNING!

Bodily injury due to tipping over of the instrument.

- ▶ Position the funnels only above the deposit spaces (see diagram).
- ▶ Maximum load per funnel holder: 2 kg.

1. Position the funnel holders above the deposit spaces.
2. Insert the funnels into the funnel holders (if necessary use the reducing inserts).
3. Position the filling container on the deposit space.
4. Fill off / filtrate the medium.

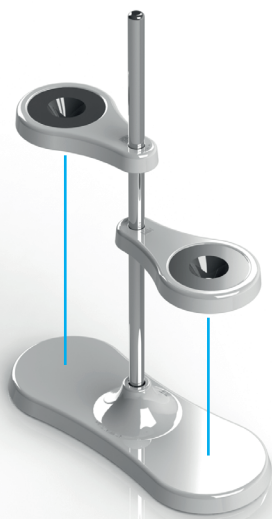Correct position of the funnel holders

Wrong position of the funnel holders -Danger of tipping over!

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