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Directions for use | Instrucciones para el uso | Instruction d'emploi | Инструкция по пользованию

Art.-Nr. 9630-0090

Pipettenständer

Pipette stand

Soporte para pipetas

Supports à pipettes

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General safety requirements

- ▶ Read these directions for use carefully before using.
- ▶ Keep these directions for use where they will be accessible for all users.
- ▶ Failure to comply with the directions for use shall render all warranty conditions null and void and may lead to personal injury and material damage.
- ▶ The instrument must only be used by qualified skilled employees.

Safety instructions before using

- ▶ Follow the manufacturer's instructions and safety instructions of the medium.
- ▶ Damage to the respiratory tract and other organs due to off-gassing media. Use only in ventilated rooms or under a suitable exhaust hood.
- ▶ The instrument must only be used within the confines of applicable material resistance and for its designated purpose. Any other type of use will affect the warranty/guarantee.
- ▶ The instrument is suitable for indoor use.
- ▶ The instrument is designed for stationary use and is not suitable for the transport of pipettes.

Safety instructions during use

WARNING!

Danger due to unsuitable working area. Danger of tipping over.

- ▶ Bright, dry and clean working place at table height without obstructions.
- ▶ Horizontal stand of the frame with sufficient space.
- ▶ The instrument has to be mounted on a horizontal, even surface.
- ▶ The device can be used in the temperature range of 0 - 40 °C.
- ▶ The instrument is suitable for the use with pipettes, thermometers and test tubes.

Layout and function

- ▶ The instrument consists of a foot, a drip pan, two perforated plates made of polypropylen (PP) and a support rod made of chrome-plated steel.
- ▶ The instrument is suitable to hold pipettes, thermometers and test tubes. It serves for storing, draining and drying pipettes.

- ▶ The instrument is suitable for 40 pipettes with a maximum diameter of 18 mm. Number and diameter of the openings:
 - 4 x Ø 18 mm
 - 4 x Ø 16 mm
 - 8 x Ø 14 mm
 - 8 x Ø 12 mm
 - 16 x Ø 10 mm
- ▶ The instrument is maintenance-free.


Scope of delivery

1 x	Support rod, Ø 16 mm, chrome-plated steel	①
1 x	Foot, PP	②
1 x	Drip pan with magnets, PP	③
2 x	Perforated plates with magnets, PP	④
1 x	Directions for use	


- ▶ Check delivery for completeness and transport damage.
- ▶ Point out any damage in the packaging to the freight forwarder/supplier immediately.

Assembly

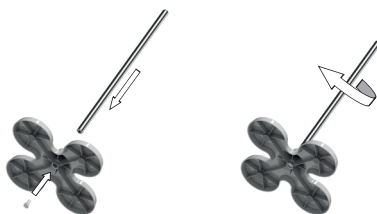1. Screw off the screw from the support rod ①.
2. Insert the support rod from above into the foot ②.
3. Fasten the support rod with the aid of the screw from below (by rotating the support rod).

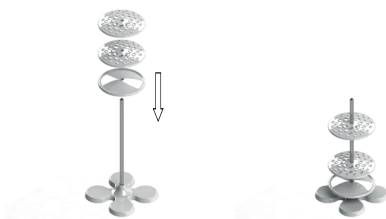4. Plug the drip pan ③ onto the support rod and slide it downwards.
5. Plug the perforated plates ④ onto the support rod and move them to the desired height.

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