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
Directions for use | Instrucciones para el uso | Instruction d'emploi | Инструкция по пользованию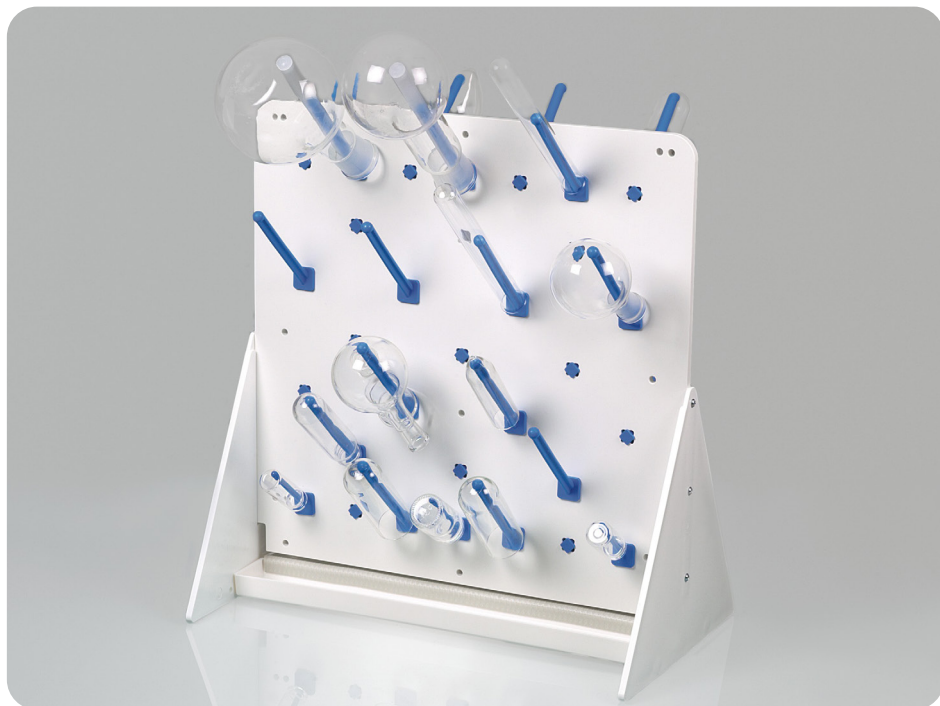

Art.-Nr. 9640-3040 / 9640-3050 / 9640-3060

Abtropfgestell

Draining rack, benchtop

Escurridor sobremesa

Egouttoir de paille

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General safety requirements

- ▶ Read these directions for use carefully before using.
- ▶ Keep these directions for use where they will be accessible for all users.
- ▶ Failure to comply with the directions for use shall render all warranty conditions null and void and may lead to personal injury and material damage.
- ▶ The instrument must only be used by qualified skilled employees.

Safety instructions before using

- ▶ The instrument must only be used within the confines of applicable material resistance and for its designated purpose. Any other type of use will affect the warranty/guarantee.
- ▶ The instrument is suitable for indoor use.
- ▶ The instrument is designed for stationary use and is not suitable for the transport of laboratory containers.
- ▶ Regularly inspect the screws and rods for tightness and integrity.

Safety instructions during use

⚠ WARNING!

Danger due to unsuitable working area. Danger of tipping over.

- ▶ Bright, dry and clean working place without obstructions.
- ▶ Horizontal stand of the frame with sufficient space to load the rack.

- ▶ The instrument has to be mounted on a horizontal, even and non-slip surface without draught.
- ▶ The device can be used in the temperature range of 0 - 60 °C.
- ▶ The instrument is not suitable for use in EX protection areas.

Layout and function

- ▶ The instrument is suitable to drain unfilled, cleaned laboratory containers with a maximum temperature of 60 °C, a one-sided maximum weight of 1.5 kg and a maximum lever arm of 30 cm.
- ▶ The instrument is maintenance-free.

Parts in contact with medium

- ▶ Draining board, connecting profile for draining channel (PVC)

- ▶ Side panel, draining channel, rods/extension rods (PP)
- ▶ Extension rod cap (LDPE)

Assembly




1. Remove the dust protection film from the surface of the board.
2. Attach the two side panels ② at the draining board ① using the provided screws ④.

CAUTION!

Risk of crushing when screwing on the side panels.

- ▶ Keep fingers away from the space between the board and the side panels.

3. Insert a bung ③ into the holes of each side panel. These serve as a stop for the draining channel ⑤.
4. Insert the rods ⑥ ⑦ into the desired holes of the draining board and lock them by turning them with the provided key ⑧ by 45°.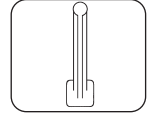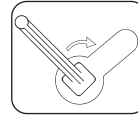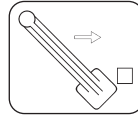

CAUTION!

Risk of crushing or scraping when mounting the rods.

- ▶ Mount the rods carefully.

CAUTION!

Bodily injury due to breakage during assembly.

- ▶ Only mount the rods, when the draining rack is not loaded.

NOTE!

The instrument is only suitable for use with original rods (see accessories).

The extension 150 mm ⑥ can only be used with rod 100 mm ⑦.

5. Push the draining channel under the draining rack up to the stop.

Safe usage

- ▶ Place one unfilled, cleaned laboratory container per rod and let it dry.
- ▶ Empty the draining channel after use.

⚠ WARNING!


Bodily injury due to tipping over of the rack or deformation of the rods.

- ▶ Maximum one-sided load per side: 1.5 kg.
- ▶ Maximum length of the laboratory containers: 30 cm.
- ▶ Maximum load per rod: 1.5 kg (rod 100 mm) and 0.8 kg (rod 60 mm).
- ▶ Pay attention to unfavorable leverage.

CAUTION!

Bodily injury due to improper use. Risk of breakage during use.

- ▶ Verify that the rods are attached correctly before use.
- ▶ Place the laboratory containers completely onto the rods.
- ▶ Choose the size of the rods according to the size of the laboratory container.
- ▶ Hang only one laboratory container per rod.
- ▶ Take special care when using laboratory glassware.

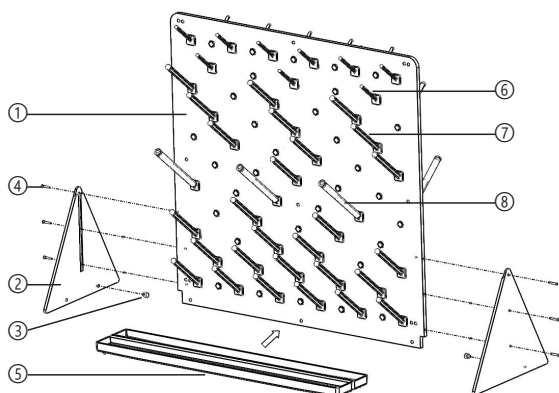**Scope of delivery**

Item no.	Board dimensions ①	Rods 60 mm ⑥	Rods 100 mm ⑦	Extension rod 150 mm ⑧	Draining Channel ⑤
9640-3040	400 x 400 mm	5	26	2	2 x 400 mm
9640-3050	500 x 500 mm	10	39	4	2 x 500 mm
9640-3060	600 x 600 mm	15	56	6	2 x 600 mm

All versions: 2 x side panel ②, 6 x screws ④, 2 x bung ③, 1 x key for rods ⑨

- ▶ Check delivery for completeness and transport damage.
- ▶ Point out any damage in the packaging to the freight forwarder/supplier immediately.

Accessories

Type	Specification	Item no.
Rods	Length 60 mm, Ø 6 mm	9640-0421
Rods	Length 100 mm, Ø 10 mm	9640-0410
Extension rods	Length 150 mm, Ø 12 mm (can only be used with 9640-0410)	9640-0415

More accessories can be found at www.buerkle.de

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