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
Directions for use | Instrucciones para el uso | Instruction d'emploi | Инструкция по пользованию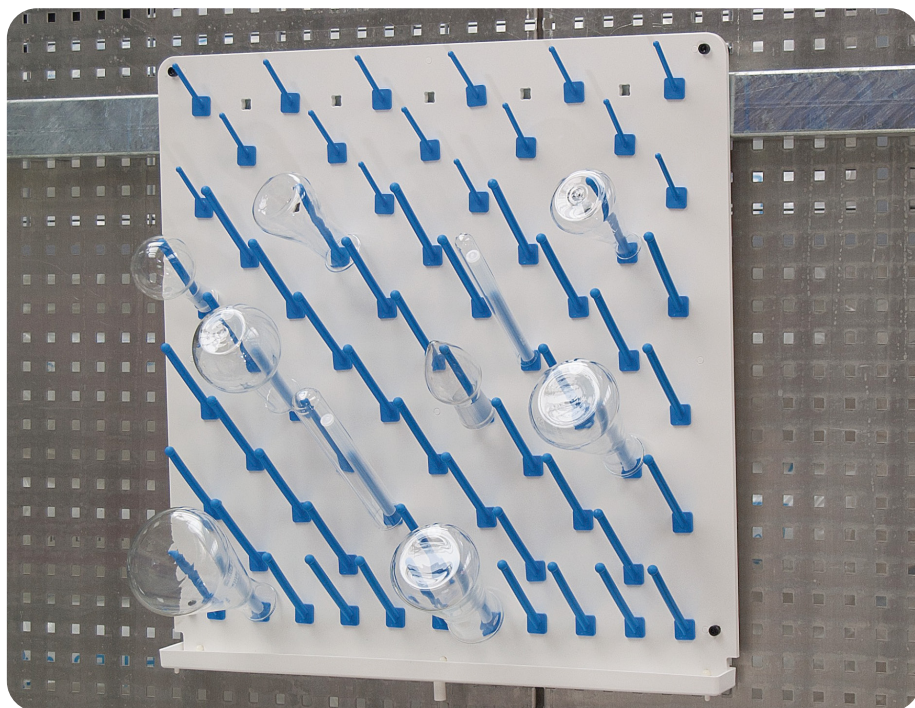

Art.-Nr. 9640-1040 / 9640-1050 / 9640-1060 / 9640-2040 / 9640-2050 / 9640-2060

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Draining rack, wallmount

Escurridor

Egouttoir mural

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General safety requirements

- ▶ Read these directions for use carefully before using.
- ▶ Keep these directions for use where they will be accessible for all users.
- ▶ Failure to comply with the directions for use shall render all warranty conditions null and void and may lead to personal injury and material damage.
- ▶ The instrument must only be used by qualified skilled employees.

Safety instructions before using

- ▶ The instrument must only be used within the confines of applicable material resistance and for its designated purpose. Any other type of use will affect the warranty/guarantee.
- ▶ The instrument is suitable for indoor use.
- ▶ The instrument is designed for stationary use and is not suitable for the transport of laboratory containers.
- ▶ Regularly inspect the screws and rods for tightness and integrity.

Safety instructions during use

⚠ WARNING!

Danger due to unsuitable working area.

- ▶ Bright, dry and clean working place without obstructions.
 - ▶ Secure suspension of the rack with sufficient space for loading.
- ▶ The rack must be hung on a suitable wall.
- ▶ The device can be used in the temperature range of 0 - 60 °C.
- ▶ The instrument is not suitable for use in EX protection areas.

Layout and function

- ▶ The instrument is suitable to drain unfilled, cleaned laboratory containers with a maximum temperature of 60 °C and a maximum lever arm of 30 cm.
- ▶ The instrument is maintenance-free.

Parts in contact with medium




- ▶ Draining board, connecting profile for draining channel (PVC)
- ▶ Side panel, draining channel, rods/extension rods (PP)
- ▶ Extension rod cap (LDPE)
- ▶ Cap nut and screw to attach the draining channel

(PA)

- ▶ Only 9640-20x0: hose (PVC), hose nozzle (LDPE)

Assembly

1. Remove the dust protection film from the surface of the board.
2. Insert the rods ① ② into the desired holes of the draining board and lock them by turning them with the provided key ④ by 45°.
3. Hold the hose nozzle ⑤ and insert the draining hose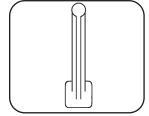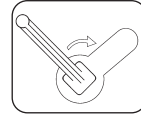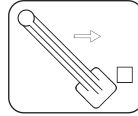

CAUTION!

Risk of crushing or scraping when mounting the rods.

- ▶ Mount the rods carefully.

CAUTION!



Bodily injury due to breakage during assembly.

- ▶ Only mount the rods, when the draining rack is not loaded.

NOTE!

The instrument is only suitable for use with original rods (see accessories).
The extension 150 mm ③ can only be used with rod 100 mm ②.

- into the nozzle (only 9640-20x0).
4. Put the draining channel ⑥ over the board and screw it on using the provided screws ⑦ and cap nuts ⑧ (only 9640-20x0).
 5. Mount the draining rack firmly to the wall using the

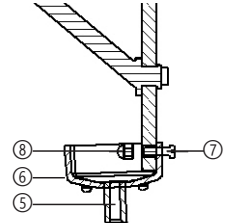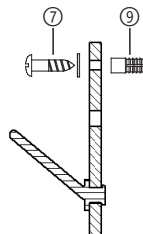provided screws ⑦ and dowels ⑨.

NOTE!

The wall mounting of the draining rack must be carried out by qualified personnel.

- ▶ The provided wall mounting set is suitable for mounting the draining rack to concrete walls or solid brick walls.
- ▶ The user himself is responsible for the correct mounting in the context of the structure of his wall.
- ▶ The minimum edge distance of the hole is 30 mm.

Safe usage

- ▶ Place one unfilled, cleaned laboratory container per rod and let it dry.
- ▶ Hang the draining hose into a laboratory basin or a suitable collecting container (only 9640-20x0).

⚠ WARNING!

Bodily injury due to deformation of the rods.

- ▶ Maximum length of the laboratory containers: 30 cm.
- ▶ Maximum load per rod: 1.5 kg (rod 100 mm) and 0.8 kg (rod 60 mm).
- ▶ Pay attention to unfavorable leverage.

WARNING!

Bodily injury due to improper use. Risk of breakage during use.

- ▶ Verify that the rods are attached correctly before use.
- ▶ Place the laboratory containers completely onto the rods.
- ▶ Choose the size of the rods according to the size of the laboratory container.
- ▶ Hang only one laboratory container per rod.
- ▶ Take special care when using laboratory glassware.

Scope of delivery

Item no.	Board dimensions	Rods 60 mm ①	Rods 100 mm ②	Extension rod 150 mm ③	Draining Channel ④
Draining rack without draining channel					
9640-1040	400 x 400 mm	5	26	2	---
9640-1050	500 x 500 mm	10	39	4	---
9640-1060	600 x 600 mm	15	56	6	---
Draining rack with draining channel					
9640-2040	400 x 400 mm	5	26	2	2 x 400 mm
9640-2050	500 x 500 mm	10	39	4	2 x 500 mm
9640-2060	600 x 600 mm	15	56	6	2 x 600 mm
All versions: 1 x wall mounting set ⑤ ⑥, 1 x key for rods ⑦					

- ▶ Check delivery for completeness and transport damage.
- ▶ Point out any damage in the packaging to the freight forwarder/supplier immediately.

Accessories

Type	Specification	Item no.
Rods	Length 60 mm , Ø 6 mm	9640-0421
Rods	Length 100 mm , Ø 10 mm	9640-0410
Extension rods	Length 150 mm , Ø 12 mm (can only be used with 9640-0410)	9640-0415

More accessories can be found at www.buerkle.de

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