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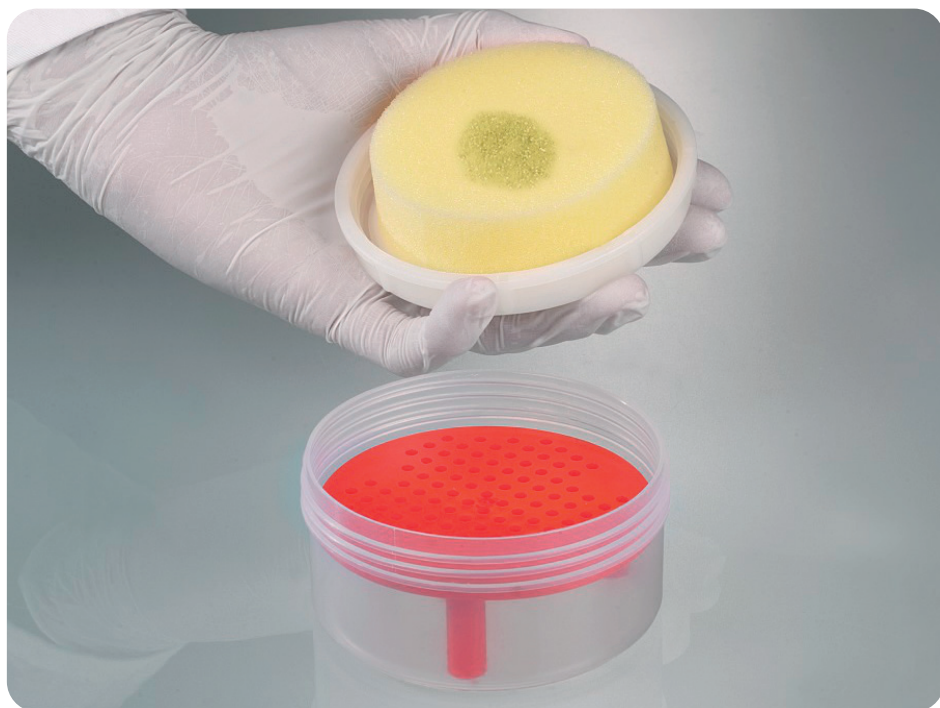
Directions for use | Instrucciones para el uso | Instruction d'emploi | Инструкция по пользованию

Art.-Nr. 9789-0001

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Mercury collector

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General safety requirements

- ▶ Read these directions for use carefully before using.
- ▶ Keep these directions for use where they will be accessible for all users.
- ▶ Failure to comply with the directions for use shall render all warranty conditions null and void and may lead to personal injury and material damage.
- ▶ The instrument must only be used by qualified skilled employees.

Safety instructions before using

- ▶ The instrument is used for safely collecting small amounts (< 100 ml) of spilled mercury.
- ▶ Follow the manufacturer's instructions and safety instructions of the chemical producer.

WARNING!

Damage to the respiratory tract and other organs due to escaping gases and liquids.

- ▶ Use the instrument only under a suitable exhaust hood and in a temperature range from 10° to 40° C.
- ▶ Wear appropriate protective clothing.

Layout and function

The mercury collector consists of a screw-on nozzle ① made of polypropylene (PP) into which a mercury separation disc ② made of PP is inserted. A special plastic foam insert ③ made of PUR polyether is used.

Point out any damage in the packaging to the freight forwarder/supplier immediately.

Safe use

CAUTION!

Before using the container, perform a visual inspection for damage and check the plastic foam insert to ensure it is properly seated and in good condition.

- ▶ If there is any damage, replace before using.
- ▶ Contact the dealer or manufacturer.

WARNING!

Damage to the respiratory tract and other organs due to escaping gases and liquids.

- ▶ Use the instrument only under a suitable exhaust hood and in a temperature range from 10° to 40° C.
- ▶ Wear appropriate protective clothing.

1. Collect the spilled mercury by lightly dabbing with the cover.
2. Press or turn the cover to wipe off the mercury on the separation disc. (Steps 1 to 2 can be repeated until all of the mercury is collected.)
3. Close the tub securely after collecting the mercury.
4. Hold the tub horizontal and set it down on a straight, non-slip base.
5. Put the mercury you collected back in the original container or dispose of it properly.

CAUTION!

The disc insert is loose. It can fall out and squirt mercury as it does so.

- ▶ Always hold the tub horizontal.

- ▶ Do not store the mercury in the tub.
- ▶ For brief transport, close the tub securely.

After usage

- ▶ Empty the mercury collector after use until there is no residue.
- ▶ When the mercury collector is placed in storage, there must be no more residue of mercury present.
- ▶ If necessary replace the plastic foam insert.

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